

## 线粒体膜电位检测试剂盒(JC-1)

产品编号	产品名称	包装
C2006	线粒体膜电位检测试剂盒(JC-1)	>100次

### 产品简介:

- 线粒体膜电位检测试剂盒(JC-1) (Mitochondrial membrane potential assay kit with JC-1)是一种以JC-1为荧光探针,快速灵敏地检测细胞、组织或纯化的线粒体膜电位变化的试剂盒,可以用于早期的细胞凋亡检测。
- JC-1是一种广泛用于检测线粒体膜电位(mitochondrial membrane potential)  $\Delta\Psi_m$ 的理想荧光探针。可以检测细胞、组织或纯化的线粒体膜电位。在线粒体膜电位较高时,JC-1聚集在线粒体的基质(matrix)中,形成聚合物(J-aggregates),可以产生红色荧光;在线粒体膜电位较低时,JC-1不能聚集在线粒体的基质中,此时JC-1为单体(monomer),可以产生绿色荧光。这样就可以非常方便地通过荧光颜色的转变来检测线粒体膜电位的变化。常用红绿荧光的相对比例来衡量线粒体去极化的比例。
- 线粒体膜电位的下降是细胞凋亡早期的一个标志性事件。通过JC-1从红色荧光到绿色荧光的转变可以很容易地检测到细胞膜电位的下降,同时也可以利用JC-1从红色荧光到绿色荧光的转变作为细胞凋亡早期的一个检测指标。JC-1检测线粒体膜电位的效果参考图1。

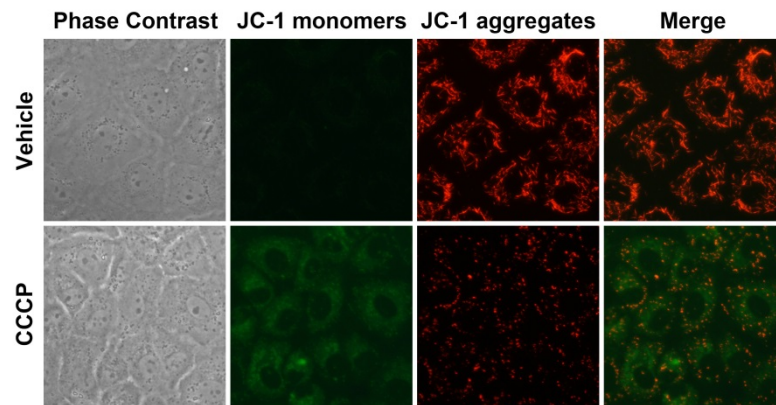


图1. 本试剂盒检测NRK-52E细胞(大鼠肾小管上皮细胞)线粒体膜电位的效果图。正常的NRK-52E细胞线粒体中JC-1以聚合物形式存在,呈明亮的红色荧光,细胞中的绿色荧光非常弱;使用CCCP处理使线粒体膜电位下降后,JC-1便不能以聚合物形式存在线粒体基质中,此时线粒体内红色荧光强度显著降低,而细胞浆中的绿色荧光显著增强。

- JC-1单体的最大激发波长为514nm,最大发射波长为529nm;JC-1聚合物(J-aggregates)的最大激发波长为585nm,最大发射波长为590nm。实际观察时,使用常规的观察红色荧光和绿色荧光的设置即可。JC-1单体和聚合物的激发光和发射光光谱参考图2。

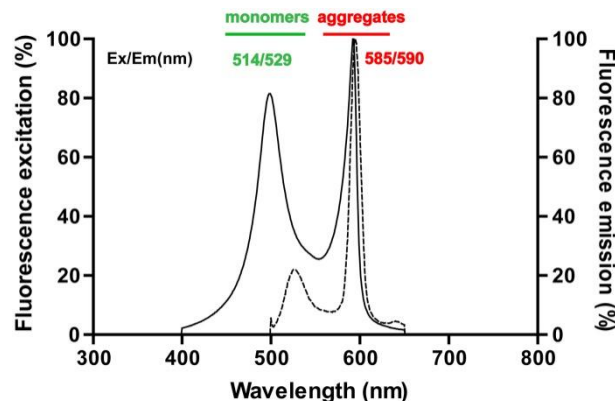


图2. JC-1单体和聚合物的激发光和发射光光谱。

- 本试剂盒提供了CCCP作为诱导线粒体膜电位下降的阳性对照。
- 对于六孔板中的样品,本试剂盒共可以检测100个样品;对于12孔中的样品,本试剂盒共可以检测200个样品。

### 包装清单:

产品编号	产品名称	包装
C2006-1	JC-1(200X)	100 $\mu$ l/管,共5管

C2006-2	超纯水	90ml
C2006-3	JC-1染色缓冲液(5X)	80ml
C2006-4	CCCP(10mM)	20 $\mu$ l
—	说明书	1份

### 保存条件：

-20°C保存。JC-1(200X)需避光保存，并尽量避免反复冻融。超纯水和JC-1染色缓冲液(5X)也可4°C保存。

### 注意事项：

- JC-1(200X)在4°C、冰浴等较低温度情况下会凝固而粘在离心管管底、管壁或管盖内，可以20-25°C水浴温育片刻至全部融解后使用。
- 必须先把JC-1(200X)用试剂盒提供的超纯水充分溶解混匀后，才可以加入JC-1染色缓冲液(5X)。不可先配制JC-1染色缓冲液(1X)再加入JC-1(200X)，这样JC-1会很难充分溶解，会严重影响后续的检测。
- 装载完JC-1后用JC-1染色缓冲液(1X)洗涤时，使JC-1染色缓冲液(1X)保持4°C左右，此时的洗涤效果较好。
- JC-1探针装载完并洗涤后尽量在30分钟内完成后续检测。在检测前需冰浴保存。
- 请勿把JC-1染色缓冲液(5X)全部配制成JC-1染色缓冲液(1X)，本试剂盒使用过程中需直接使用JC-1染色缓冲液(5X)。
- 如果发现JC-1染色缓冲液(5X)中有沉淀，必须全部溶解后才能使用，为促进溶解可以在37°C加热。
- CCCP为线粒体电子传递链抑制剂，对人体有害，操作时请小心，并注意有效防护以避免直接接触人体或吸入体内。
- 本产品仅限于专业人员的科学研究用，不得用于临床诊断或治疗，不得用于食品或药品，不得存放于普通住宅内。
- 为了您的安全和健康，请穿实验服并戴一次性手套操作。

### 使用说明：

#### 1. JC-1染色工作液的配制：

六孔板每孔所需JC-1染色工作液的量为1ml，其它培养器皿的JC-1染色工作液的用量以此类推；对于细胞悬液每50-100万细胞需0.5ml JC-1染色工作液。取适量JC-1(200X)，按照每50 $\mu$ l JC-1(200X)加入8ml超纯水的比例稀释JC-1。剧烈Vortex充分溶解并混匀JC-1。然后再加入2ml JC-1染色缓冲液(5X)，混匀后即为JC-1染色工作液。

#### 2. 阳性对照的设置：

把试剂盒中提供的CCCP(10mM)推荐按照1:1000的比例加入到细胞培养液中，稀释至10 $\mu$ M，处理细胞20分钟。随后按照下述方法装载JC-1，进行线粒体膜电位的检测。对于大多数细胞，通常10 $\mu$ M CCCP处理20分钟后线粒体的膜电位会完全丧失，JC-1染色后观察应呈绿色荧光；而正常的细胞经JC-1染色后应显示红色荧光。对于特定的细胞，CCCP的作用浓度和作用时间可能有所不同，需自行参考相关文献资料确定。

#### 3. 对于悬浮细胞：

- a. 取10-60万细胞，重悬于0.5ml细胞培养液中，细胞培养液中可以含血清和酚红。
- b. 加入0.5ml JC-1染色工作液，颠倒数次混匀。细胞培养箱中37°C孵育20分钟。
- c. 在孵育期间，按照每1ml JC-1染色缓冲液(5X)加入4ml蒸馏水的比例，配制适量的JC-1染色缓冲液(1X)，并放置于冰浴。
- d. 37°C孵育结束后，600g 4°C离心3-4分钟，沉淀细胞。弃上清，注意尽量不要吸除细胞。
- e. 用JC-1染色缓冲液(1X)洗涤2次：加入1ml JC-1染色缓冲液(1X)重悬细胞，600g 4°C离心3-4分钟，沉淀细胞，弃上清。再加入1ml JC-1染色缓冲液(1X)重悬细胞，600g 4°C离心3-4分钟，沉淀细胞，弃上清。
- f. 再用适量JC-1染色缓冲液(1X)重悬后，用荧光显微镜或激光共聚焦显微镜观察，也可以用荧光分光光度计检测或流式细胞仪分析。

#### 4. 对于贴壁细胞：

**注意：**对于贴壁细胞，如果希望采用荧光分光光度计或流式细胞仪检测，可以先收集细胞，重悬后参考悬浮细胞的检测方法。

- a. 对于六孔板的一个孔，吸除培养液，根据具体实验如有必要可以用PBS或其它适当溶液洗涤细胞一次，加入1ml细胞培养液。细胞培养液中可以含有血清和酚红。
- b. 加入1ml JC-1染色工作液，充分混匀。细胞培养箱中37°C孵育20分钟。
- c. 在孵育期间，按照每1ml JC-1染色缓冲液(5X)加入4ml蒸馏水的比例，配制适量的JC-1染色缓冲液(1X)，并放置于冰浴。
- d. 37°C孵育结束后，吸除上清，用JC-1染色缓冲液(1X)洗涤2次。
- e. 加入2ml细胞培养液，培养液中可以含有血清和酚红。
- f. 荧光显微镜或激光共聚焦显微镜下观察。

#### 5. 对于纯化的线粒体：

- a. 把配制好的JC-1染色工作液再用JC-1染色缓冲液(1X)稀释5倍。
- b. 0.9ml 5倍稀释的JC-1染色工作液中加入0.1ml总蛋白量为10-100 $\mu$ g纯化的线粒体。
- c. 用荧光分光光度计或荧光酶标仪检测：混匀后直接用荧光分光光度计进行时间扫描(time scan)，激发波长为485nm，发射波长为590nm。如果使用荧光酶标仪，激发波长不能设置为485nm时，可以在475-520nm范围内设置激发波长。另外，也可以参考下面步骤6中的波长设置进行荧光检测。
- d. 用荧光显微镜或激光共聚焦显微镜观察：方法同下面的步骤6。

#### 6. 荧光观测和结果分析：

- a. 检测JC-1单体时可以把激发光设置为490nm，发射光设置为530nm；检测JC-1聚合物时，可以把激发光设置为525nm，发射光设置为590nm。注意：此处测定荧光时不必把激发光和发射光设置在最大激发波长和最大发射波长。如使用荧光显微镜观察，检测JC-1单体时可以参考观察其它绿色荧光时的设置，如观察GFP或FITC时的设置；检测JC-1聚合物时可以参考观察其它红色荧光，如碘化丙啶或Cy3时的设置。出现绿色荧光说明线粒体膜电位下降，并且该细胞很可能处于细胞凋亡早期。出现红色荧光说明线粒体膜电位比较正常，细胞的状态也比较正常。

#### 相关产品：

产品编号	产品名称	包装
C1002	DAPI	5mg/ml×0.2ml
C1005/C1006	DAPI 染色液	10ml/50ml
C1011	Hoechst 33258	10mg
C1017/C1018	Hoechst 33258 染色液	10ml/50ml
C1022	Hoechst 33342	10mg
C1025/C1026	Hoechst 33342 染色液	10ml/50ml
C1027/C1028	Hoechst 33342 活细胞染色液(100X)	0.1ml/0.5ml
C1031	CFDA SE (细胞增殖示踪荧光探针)	5mg
C1033	Actin-Tracker Green (微丝绿色荧光探针)	0.2ml
C1036	DiI (细胞膜红色荧光探针)	10mg
C1038	DiO (细胞膜绿色荧光探针)	10mg
C1041	ER-Tracker Red (内质网红色荧光探针)	20μl
C1043	Golgi-Tracker Red (高尔基体红色荧光探针)	1mg
C1046	Lyso-Tracker Red (溶酶体红色荧光探针)	50μl
C1048	Mito-Tracker Green (线粒体绿色荧光探针)	50μg
C1050	Tubulin-Tracker Red (微管红色荧光探针)	40μl
C2005	JC-1	1mg
C2006	线粒体膜电位检测试剂盒(JC-1)	>100 次
C2007	Rhodamine 123	5mg
S0019	DAF-FM DA (NO 荧光探针)	>100 次
S0063	Dihydroethidium (超氧化物阴离子荧光探针)	5mg
S1006	BCECF AM (pH 荧光探针, 5mM)	50 微升
S1052	Fura-2 AM (钙离子荧光探针, 2mM)	50 微升
S1056	Fluo-3 AM (钙离子荧光探针, 5mM)	20 微升
S1060	Fluo-4 AM (钙离子荧光探针, 2mM)	25 微升
S1082	MQAE (氯离子荧光探针)	20mg

#### 使用本产品的文献：

- Su X, Zheng X, Ni J. Lanthanum citrate induces anoikis of HeLa cells. *Cancer Lett.* 2009 Nov 28;285(2):200-9.
- Zhu XJ, Shi Y, Peng J, Guo CS, Shan NN, Qin P, Ji XB, Hou M. The effects of BAFF and BAFF-R-Fc fusion protein in immune thrombocytopenia. *Blood.* 2009 Dec 17;114(26):5362-7.
- Xu HN, Huang WD, Cai Y, Ding M, Gu JF, Wei N, Sun LY, Cao X, Li HG, Zhang KJ, Liu XR, Liu XY. HCCS1-armed, quadruple-regulated oncolytic adenovirus specific for liver cancer as a cancer targeting gene-viro-therapy strategy. *Mol Cancer.* 2011 Nov 1;10:133.
- Wei A, Zhou D, Xiong C, Cai Y, Ruan J. A novel non-aromatic B-ring flavonoid: isolation, structure elucidation and its induction of apoptosis in human colon HT-29 tumor cell via the reactive oxygen species-mitochondrial dysfunction and MAPK activation. *Food Chem Toxicol.* 2011 Sep;49(9):2445-52.
- Nie Y, Han BM, Liu XB, Yang JJ, Wang F, Cong XF, Chen X. Identification of MicroRNAs involved in hypoxia- and serum deprivation-induced apoptosis in mesenchymal stem cells. *Int J Biol Sci.* 2011;7(6):762-8.
- Gao M, Wang J, Wang W, Liu J, Wong CW. Phosphatidylinositol 3-kinase affects mitochondrial function in part through inducing peroxisome proliferator-activated receptor  $\gamma$  coactivator-1 $\beta$  expression. *BRIT J PHARMACOL.* 2011 Feb;162(4):1000-8.
- Liu J, Xu X, Feng X, Zhang B, Wang J. Adenovirus-mediated delivery of bFGF small interfering RNA reduces STAT3 phosphorylation and induces the depolarization of mitochondria and apoptosis in glioma cells U251. *J EXP CLIN CANC RES.* 2011 Sep 9;30:80.
- Liu Q, Li Y, Hu L, Wang D. Lycium barbarum Polysaccharides Attenuate Cisplatin-Induced Hair Cell Loss in Rat Cochlear Organotypic Cultures. *Int J Mol Sci.* 2011;12(12):8982-92.
- Gu L, Liang X, Wang L, Yan Y, Ni Z, Dai H, Gao J, Mou S, Wang Q, Chen X, Wang L, Qian J. Functional metabotropic glutamate receptors 1 and 5 are expressed in murine podocytes. *Kidney Int.* 2012 Mar;81(5):458-68.
- Ma J, Zhang L, Han W, Shen T, Ma C, Liu Y, Nie X, Liu M, Ran Y, Zhu D. Activation of JNK/c-Jun is required for the proliferation, survival, and angiogenesis induced by EET in pulmonary artery endothelial cells. *J Lipid Res.* 2012 Jun;53(6):1093-105.
- Ou XH, Li S, Wang ZB, Li M, Quan S, Xing F, Guo L, Chao SB, Chen Z, Liang XW, Hou Y, Schatten H, Sun QY. Maternal insulin resistance causes oxidative stress and mitochondrial dysfunction in mouse oocytes. *Hum Reprod.* 2012 Jul;27(7):2130-45.
- Tao Q, Fan X, Li T, Tang Y, Yang D, Le W. Gender segregation in gene expression and vulnerability to oxidative stress induced injury in ventral mesencephalic cultures of dopamine neurons. *J Neurosci Res.* 2012 Jan;90(1):167-78.

13. Shi ZY, Li YQ, Kang YH, Hu GQ, Huang-fu CS, Deng JB, Liu B. Piperonal ciprofloxacin hydrazone induces growth arrest and apoptosis of humanhepatocarcinoma SMMC-7721 cells. *Acta Pharmacol Sin.* 2012 Feb;33(2):271-8.
14. Sang H, Zhang L, Li J. Anti-benzopyrene-7,8-diol-9,10-epoxide induces apoptosis via mitochondrial pathway in humanbronchiolar epithelium cells independent of the mitochondria permeability transition pore. *Food Chem Toxicol.* 2012 Jul;50(7):2417-23.
15. Fu J, Dang Z, Deng Y, Lu G. Regulation of c-Myc and Bcl-2 induced apoptosis of human bronchial epithelial cells by zinc oxide nanoparticles. *J BIOMED NANOTECHNOL.* 2012 Aug;8(4):669-75.
16. Zhang J, Song W, Guo J, Zhang J, Sun Z, Ding F, Gao M. Toxic effect of different ZnO particles on mouse alveolar macrophages. *J HAZARD MATER.* 2012 Jun 15;219-220:148-55.
17. Khan M, Ding C, Rasul A, Yi F, Li T, Gao H, Gao R, Zhong L, Zhang K, Fang X, Ma T. Isoalantolactone induces reactive oxygen species mediated apoptosis in pancreatic carcinomaPANC-1 cells. *Int J Biol Sci.* 2012;8(4):533-47.
18. Ding F, Shao ZW, Yang SH, Wu Q, Gao F, Xiong LM. Role of mitochondrial pathway in compression-induced apoptosis of nucleus pulposus cells. *Apoptosis.* 2012 Jun;17(6):579-90.
19. Lu YS, Huang BT, Huang YX. Reactive oxygen species formation and apoptosis in human peripheral blood mononuclear cell induced by 900MHz mobile phone radiation. *Oxid Med Cell Longev.* 2012;2012:740280.
20. Li XY, Jing CQ, Lei WL, Li J, Wang JJ. Apoptosis caused by imidazolium-based ionic liquids in PC12 cells. *ECOTOX ENVIRON SAFE.* 2012 Sep;83:102-7.
21. Mo R, Sun Q, Li N, Zhang C. Intracellular delivery and antitumor effects of pH-sensitive liposomes based on zwitterionic oligopeptide lipids. *Biomaterials.* 2013 Apr;34(11):2773-86.
22. Zhang F, Kong DS, Zhang ZL, Lei N, Zhu XJ, Zhang XP, Chen L, Lu Y, Zheng SZ. Tetramethylpyrazine induces G0/G1 cell cycle arrest and stimulates mitochondrial-mediated and caspase-dependent apoptosis through modulating ERK/p53 signaling in hepatic stellate cells in vitro. *Apoptosis.* 2013 Feb;18(2):135-49.
23. Chen G, Zhang P, Huang T, Yu W, Lin J, Li P, Chen K. Polysaccharides from *Rhizopus nigricans* mycelia induced apoptosis and G2/M arrest in BGC-823 cells. *Carbohydrate Polymers.* 2013 Sep;97(2):800–8.
24. Sun J, Bai S, Bai W, Zou F, Zhang L, Li G, Hu Y, Li M, Yan R, Su Z, Huang Y. 1,3-Dichloro-2-propanol inhibits progesterone production through the expression of steroidogenic enzymes and cAMP concentration in Leydig cells. *Food Chem.* 2014 Jul 1;154:330-6.
25. Yang L, Di G, Qi X, Qu M, Wang Y, Duan H, Danielson P, Xie L, Zhou Q. Substance P promotes diabetic corneal epithelial wound healing through molecular mechanisms mediated via the neurokinin-1 receptor. *Diabetes.* 2014 Dec;63(12):4262-74.
26. Chu JH, Zhao CR, Song ZY, Wang RQ, Qin YZ, Li WB, Qu XJ. 1082-39, an analogue of sorafenib, inhibited human cancer cell growth more potently than sorafenib. *Biomed Pharmacother.* 2014 Apr;68(3):335-41.
27. Wang W, Xie Q, Zhou X, Yao J, Zhu X, Huang P, Zhang L, Wei J, Xie H, Zhou L, Zheng S. Mitofusin-2 triggers mitochondria Ca<sup>2+</sup> influx from the endoplasmic reticulum to induce apoptosis in hepatocellular carcinoma cells. *Cancer Lett.* 2015 Mar 1;358(1):47-58.
28. Zong Y, Huang Y, Chen S, Zhu M, Chen Q, Feng S, Sun Y, Zhang Q, Tang C, Du J, Jin H. Downregulation of Endogenous Hydrogen Sulfide Pathway Is Involved in Mitochondrion-Related Endothelial Cell Apoptosis Induced by High Salt. *Oxid Med Cell Longev.* 2015;2015:754670.
29. Zhang J, Wu D, Xing Z, Liang S, Han H, Shi H, Zhang Y, Yang Y, Li Q. N-Isopropylacrylamide-modified polyethylenimine-mediated p53 gene delivery to prevent the proliferation of cancer cells. *COLLOID SURFACE B.* 2015 May 1;129:54-62.
30. Wang AT, Liang DS, Liu YJ, Qi XR. Roles of ligand and TPGS of micelles in regulating internalization, penetration and accumulation against sensitive or resistant tumor and therapy for multidrug resistant tumors. *Biomaterials.* 2015 Jun;53:160-72.
31. Bai L, Guo Y, Du Y, Wang H, Zhao Z, Huang Y, Tang J. 47kDa isoform of Annexin A7 affecting the apoptosis of mouse hepatocarcinoma cells line. *Biomed Pharmacother.* 2016 Oct;83:1127-1131.
32. Zhang F, Zhu X, Gong J, Sun Y, Chen D, Wang J, Wang Y, Guo M, Li W. Lysosome-mitochondria-mediated apoptosis specifically evoked in cancer cells induced by goldnanorods. *NANOMEDICINE-UK.* 2016 Aug;11(15):1993-2006.
33. Dong K, Yan Y, Wang P, Shi X, Zhang L, Wang K, Xing J, Dong Y. Biodegradable mixed MPEG-SS-2SA/TPGS micelles for triggered intracellular release of paclitaxel andreversing multidrug resistance. *INT J NANOMED.* 2016 Oct 6;11:5109-5123.
34. Wang M, Li Y, HuangFu M, Xiao Y, Zhang T, Han M, Xu D, Li F, Ling D, Jin Y, Gao J. Pluronic-attached polyamidoamine dendrimer conjugates overcome drug resistance in breast cancer. *NANOMEDICINE-UK.* 2016 Nov;11(22):2917-2934.
35. Zhang A, Sheng Y, Zou M. Antiproliferative activity of Alisol B in MDA-MB-231 cells is mediated by apoptosis, dysregulation ofmitochondrial functions, cell cycle arrest and generation of reactive oxygen species. *Biomed Pharmacother.* 2017 Mar;87:110-117.
36. Su S, Zhu X, Lin L, Chen X, Wang Y, Zi J, Dong Y, Xie Y, Zhu Y, Zhang J, Zhu J, Xu D, Xu N, Lou X, Liu S. Lowering Endogenous Cathepsin D Abundance Results in Reactive Oxygen SpeciesAccumulation and Cell Senescence. *Mol Cell Proteomics.* 2017 Jul;16(7):1217-1232.
37. Xu D, Chen L, Chen X, Wen Y, Yu C, Yao J, Wu H, Wang X, Xia Q, Kong X. The triterpenoid CDDO-imidazolide ameliorates mouse liver ischemia-reperfusion injury through activating the Nrf2/HO-1 pathway enhanced autophagy. *Cell Death Dis.* 2017 Aug 10;8(8):e2983.
38. Huang HL, Shi YP, He HJ, Wang YH, Chen T, Yang LW, Yang T, Chen J, Cao J, Yao WM, Liu G. MiR-4673 Modulates Paclitaxel-Induced Oxidative Stress and Loss of Mitochondrial MembranePotential by Targeting 8-Oxoguanine-DNA Glycosylase-1. *CELL PHYSIOL BIOCHEM.* 2017;42(3):889-900.
39. Zhang A, Sheng Y, Zou M. Antiproliferative activity of Alisol B in MDA-MB-231 cells is mediated by apoptosis, dysregulation of mitochondrial functions, cell cycle arrest and generation of reactive oxygen species. *Biomed Pharmacother.* 2017 Mar;87:110-117.
40. Wang T, Gao YY, Chen L, Nie ZW, Cheng W, Liu X, Schatten H4, Zhang X, Miao YL. Melatonin prevents postovulatory oocyte aging and promotes subsequent embryonicdevelopment in the pig. *AGING-US.* 2017 Jun 26;9(6):1552-1564.
41. You P, Wu H, Deng M, Peng J, Li F, Yang Y. Brevilin A induces apoptosis and autophagy of colon adenocarcinoma cell CT26 viamitochondrial pathway and PI3K/AKT/mTOR inactivation. *Biomed Pharmacother.* 2017 Dec 28;98:619-625.
42. Yang X, Tang S, Li D, Yu X, Wang F, Xiao X. DIDS inhibits overexpression BAK1-induced mitochondrial apoptosis throughGSK3 β / β -catenin signaling pathway. *J Cell Physiol.* 2017 Dec 12.
43. Wang P, Gao YM, Sun X, Guo N, Li J, Wang W, Yao LP, Fu YJ. Hepatoprotective effect of 2'-O-galloylhyperin against oxidative stress-induced liver damage through induction of Nrf2/ARE-mediated antioxidant pathway. *Food Chem Toxicol.* 2017 Apr;102:129-142.
44. Shan M, Qin J, Jin F, Han X, Guan H, Li X, Zhang J, Zhang H, Wang Y. Autophagy suppresses isoprenaline-induced M2 macrophage polarization via the ROS/ERK and mTOR signaling pathway. *FREE RADICAL BIO MED.* 2017 Sep;110:432-443.
45. Huo X, Wang C, Yu Z, Peng Y, Wang S, Feng S, Zhang S, Tian X, Sun C, Liu K, Deng S, Ma X. Human transporters, PEPT1/2, facilitate melatonin transportation into mitochondria of cancercells: An implication of the therapeutic potential. *J Pineal Res.* 2017 May;62(4).
46. Yan Y, Wang L, He J, Liu P, Lv X, Zhang Y, Xu X, Zhang L, Zhang Y. Synergy with interferon-lambda 3 and sorafenib suppresses hepatocellular carcinomaproiferation. *Biomed Pharmacother.* 2017 Apr;88:395-402.
47. Zhang Q, Wu D, Yang Y, Liu T, Liu H. Dexmedetomidine Alleviates Hyperoxia-Induced Acute Lung Injury via Inhibiting NLRP3Inflammasome

- Activation. *CELL PHYSIOL BIOCHEM*. 2017;42(5):1907-1919.
48. Yang J, Yu J, Li D, Yu S, Ke J, Wang L, Wang Y, Qiu Y, Gao X, Zhang J, Huang L. Store-operated calcium entry-activated autophagy protects EPC proliferation via the CAMKK2-MTOR pathway in ox-LDL exposure. *Autophagy*. 2017 Jan 2;13(1):82-98.
  49. Hu F, Zhao Y, Yu Y, Fang JM, Cui R, Liu ZQ, Guo XL, Xu Q. Docetaxel-mediated autophagy promotes chemoresistance in castration-resistant prostate cancer cells by inhibiting STAT3. *Cancer Lett*. 2018 Mar 1;416:24-30.
  50. Jiang HY, Yang Y, Zhang YY, Xie Z, Zhao XY, Sun Y, Kong WJ. The dual role of poly(ADP-ribose) polymerase-1 in modulating parthanatos and autophagy under oxidative stress in rat cochlear marginal cells of the stria vascularis. *Redox Biol*. 2018 Apr;14:361-370.
  51. Wang Z, Zhou F, Dou Y, Tian X, Liu C, Li H, Shen H, Chen G. Melatonin Alleviates Intracerebral Hemorrhage-Induced Secondary Brain Injury in Rats via Suppressing Apoptosis, Inflammation, Oxidative Stress, DNA Damage, and Mitochondria Injury. *Transl Stroke Res*. 2018 Feb;9(1):74-91.
  52. Zhang H, Chen S, Zeng M, Lin D, Wang Y, Wen X, Xu C, Yang L, Fan X, Gong Y, Zhang H, Kong X. Apelin-13 Administration Protects Against LPS-Induced Acute Lung Injury by Inhibiting NF- $\kappa$ B Pathway and NLRP3 Inflammasome Activation. *CELL PHYSIOL BIOCHEM*. 2018;49(5):1918-1932.
  53. Xue P, Yang R, Sun L, Li Q, Zhang L, Xu Z, Kang Y. Indocyanine Green-Conjugated Magnetic Prussian Blue Nanoparticles for Synchronous Photothermal/Photodynamic Tumor Therapy. *NANO-MICRO LETT*. 2018;10(4):74.
  54. Li S, Yang X, Feng Z, Wang P, Zhu W, Cui S. Catalase Enhances Viability of Human Chondrocytes in Culture by Reducing Reactive Oxygen Species and Counteracting Tumor Necrosis Factor- $\alpha$ -Induced Apoptosis. *CELL PHYSIOL BIOCHEM*. 2018;49(6):2427-2442.
  55. Hu Y, Yu K, Wang G, Zhang D, Shi C, Ding Y, Hong D, Zhang D, He H, Sun L, Zheng JN, Sun S, Qian F. Lanatoside C inhibits cell proliferation and induces apoptosis through attenuating Wnt/ $\beta$ -catenin/c-Myc signaling pathway in human gastric cancer cell. *Biochem Pharmacol*. 2018 Apr;150:280-292.
  56. Jiang HY, Yang Y, Zhang YY, Xie Z, Zhao XY, Sun Y, Kong WJ. The dual role of poly(ADP-ribose) polymerase-1 in modulating parthanatos and autophagy under oxidative stress in rat cochlear marginal cells of the stria vascularis. *Redox Biol*. 2018 Apr;14:361-370.
  57. Liu X, Zhang N, Wang D, Zhu D, Yuan Q, Zhang X, Qian L, Niu H, Lu Y, Ren G, Tian K, Yuan H. Downregulation of reticulocalbin-1 differentially facilitates apoptosis and necroptosis in human prostate cancer cells. *Cancer Sci*. 2018 Apr;109(4):1147-1157.
  58. Peng J, Zhou Y, Zhang Z, Wang Z, Gao L, Zhang X, Fang Z, Li G, Chen H, Yang H, Gao L. The detrimental effects of glucocorticoids exposure during pregnancy on offspring's cardiac functions mediated by hypermethylation of bone morphogenetic protein-4. *Cell Death Dis*. 2018 Aug 6;9(8):834.
  59. Gong Sun X, Zhao Y, Jiang B, Xin Z, Shi M, Song L, Qin Q, Wang Q, Liu X. Inhibition of MUC1-C regulates metabolism by AKT pathway in esophageal squamous cell carcinoma. *J Cell Physiol*. 2018 Dec 6.
  60. You P, Wu H, Deng M, Peng J, Li F, Yang Y. Brevilin A induces apoptosis and autophagy of colon adenocarcinoma cell CT26 via mitochondrial pathway and PI3K/AKT/mTOR inactivation. *Biomed Pharmacother*. 2018 Feb;98:619-625.
  61. Wang Z, Zhou F, Dou Y, Tian X, Liu C, Li H, Shen H, Chen G. Melatonin Alleviates Intracerebral Hemorrhage-Induced Secondary Brain Injury in Rats via Suppressing Apoptosis, Inflammation, Oxidative Stress, DNA Damage, and Mitochondria Injury. *Transl Stroke Res*. 2018 Feb;9(1):74-91.
  62. Maryam A, Mehmood T, Yan Q, Li Y, Khan M, Ma T. Proscillaridin A Promotes Oxidative Stress and ER Stress, Inhibits STAT3 Activation, and Induces Apoptosis in A549 Lung Adenocarcinoma Cells. *Oxid Med Cell Longev*. 2018 Jan 11;2018:3853409.
  63. Qi G, Guo R, Tian H, Li L, Liu H, Mi Y, Liu X. Nobiletin protects against insulin resistance and disorders of lipid metabolism by reprogramming of circadian clock in hepatocytes. *BBA-MOL CELL BIOL L*. 2018 Jun;1863(6):549-562.
  64. Yang X, Tang S, Li D, Yu X, Wang F, Xiao X. DIDS inhibits overexpression BAK1-induced mitochondrial apoptosis through GSK3 $\beta$ / $\beta$ -catenin signaling pathway. *J Cell Physiol*. 2018 Jun;233(6):5070-5077.
  65. Han H, Chen W, Yang J, Liang X, Wang Y, Li Q, Yang Y, Li K. Inhibition of cell proliferation and migration through nucleobase-modified polyamidoamine-mediated p53 delivery. *INT J NANOMED*. 2018 Mar 6;13:1297-1311.
  66. Zhao Y, Zhao R, Wu J, Wang Q, Pang K, Shi Q, Gao Q, Hu Y, Dong X, Zhang J, Sun J. Melatonin protects against A $\beta$ -induced neurotoxicity in primary neurons via miR-132/PTEN/AKT/FOXO3a pathway. *Biofactors*. 2018 Nov;44(6):609-618.
  67. Wu S, Lei L, Song Y, Liu M, Lu S, Lou D, Shi Y, Wang Z, He D. Mutation of hop-1 and pink-1 attenuates vulnerability of neurotoxicity in *C. elegans*: the role of mitochondria-associated membrane proteins in Parkinsonism. *Exp Neurol*. 2018 Nov;309:67-78.
  68. Liu HN, Guo NN, Guo WW, Huang-Fu MY, Vakili MR, Chen JJ, Xu WH, Wei QC, Han M, Lavasanifar A, Gao JQ. Delivery of mitochondriotropic doxorubicin derivatives using self-assembling hyaluronic acid nanocarriers in doxorubicin-resistant breast cancer. *Acta Pharmacol Sin*. 2018 Oct;39(10):1681-1692.
  69. Liu L, Zuo Z, Lu S, Wang L, Liu A, Liu X. Silencing of PINK1 represses cell growth, migration and induces apoptosis of lung cancer cells. *Biomed Pharmacother*. 2018 Oct;106:333-341.
  70. Jin L, Cai Q, Wang S, Wang S, Mondal T, Wang J, Quan Z. Long noncoding RNA MEG3 regulates LATS2 by promoting the ubiquitination of EZH2 and inhibits proliferation and invasion in gallbladder cancer. *Cell Death Dis*. 2018 Oct 3;9(10):1017.
  71. Zhang CC, Li CG, Wang YF, Xu LH, He XH, Zeng QZ, Zeng CY, Mai FY, Hu B, Ouyang DY. Chemotherapeutic paclitaxel and cisplatin differentially induce pyroptosis in A549 lung cancer cells via caspase-3/GSDME activation. *Apoptosis*. 24(3-4):312-325.
  72. Xu Z, Zhang Y, Zhang S, Jia X, Zhong G, Yang Y, Du Q, Li J, Liu Z. Novel half-sandwich iridium O<sup>+</sup>C (carbene)-Complexes: In vitro and in vivo tumor growth suppression and pro-apoptosis via ROS-mediated cross-talk between mitochondria and lysosomes. *Cancer Lett*. 447:75-85.
  73. Du X, Zhang T, Ma G, Gu X, Wang G, Li J. Glucose-responsive mesoporous silica nanoparticles to generation of hydrogen peroxide for synergistic cancer starvation and chemotherapy. *INT J NANOMED*. 14:2233-2251.
  74. Zhao SY, Liao LX, Tu PF, Li WW, Zeng KW. Icaritin Inhibits AGE-Induced Injury in PC12 Cells by Directly Targeting Apoptosis Regulator Bax. *Oxid Med Cell Longev*. 2019:7940808.
  75. Shan C, Gong YL, Zhuang QQ, Hou YF, Wang SM, Zhu Q, Huang GR, Tao B, Sun LH, Zhao HY, Li ST, Liu JM. Protective effects of  $\beta$ -nicotinamide adenine dinucleotide against motor deficits and dopaminergic neuronal damage in a mouse model of Parkinson's disease. *PROG NEURO-PSYCHOPH*. 94:109670.
  76. Wang HW, Zhang Y, Tan PP, Jia LS, Chen Y, Zhou BH. Mitochondrial respiratory chain dysfunction mediated by ROS is a primary point of fluoride-induced damage in Hepa1-6 cells. *Environ Pollut*. 255(Pt 3):113359.
  77. Song H, Wang M, Xin T. Mst1 contributes to nasal epithelium inflammation via augmenting oxidative stress and mitochondrial dysfunction in a manner dependent on Nrf2 inhibition. *J Cell Physiol*. 234(12):23774-23784.
  78. Wu M, Chen W, Zhang S, Huang S, Zhang A, Zhang Y, Jia Z. Rotenone protects against  $\beta$ -cell apoptosis and attenuates type 1 diabetes mellitus. *Apoptosis*. 24(11-12):879-891.
  79. Zhou J, Xue Z, He HN, Liu X, Yin SY, Wu DY, Zhang X, Schatten H, Miao YL. Resveratrol delays postovulatory aging of mouse oocytes through activating mitophagy. *AGING-US*. 11(23):11504-11519.
  80. Huang X, Ni B, Xi Y, Chu X, Zhang R, You H. Protease-activated receptor 2 (PAR-2) antagonist AZ3451 as a novel therapeutic agent for osteoarthritis. *AGING-US*. 11(24):12532-12545.
  81. Tong Y, Feng A, Hou X, Zhou Q, Hu X. Nanoholes Regulate the Phytotoxicity of

- Single-Layer Molybdenum Disulfide. *Environ Sci Technol*. 53(23):13938-13948.
82. Zhang S, Liu C, Zhang X. Mitochondrial Damage Mediated by miR-1 Overexpression in Cancer Stem Cells. *MOL THER-NUCL ACIDS*. 18:938-953.
83. Wang L, Yan W, Wang J. Silencing MEKK3 attenuates cardiomyocyte injury caused by hypoxia/reoxygenation via the sonic hedgehog pathway. *J Cell Physiol*.
84. Li Z, Li Q, Lv W, Jiang L, Geng C, Yao X, Shi X, Liu Y, Cao J. The interaction of Atg4B and Bcl-2 plays an important role in Cd-induced crosstalk between apoptosis and autophagy through disassociation of Bcl-2-Beclin1 in A549 cells. *FREE RADICAL BIO MED*. 130:576-591.
85. Yao X, Zhang J, Jing X, Ye Y, Guo J, Sun K, Guo F. Fibroblast growth factor 18 exerts anti-osteoarthritic effects through PI3K-AKT signaling and mitochondrial fusion and fission. *Pharmacol Res*. 139:314-324.
86. Fan Y, Yang Q, Yang Y, Gao Z, Ma Y, Zhang L, Liang W, Ding G. Sirt6 Suppresses High Glucose-Induced Mitochondrial Dysfunction and Apoptosis in Podocytes through AMPK Activation. *Int J Biol Sci*. 15(3):701-713.
87. Cai G, Yu W, Song D, Zhang W, Guo J, Zhu J, Ren Y, Kong L. Discovery of fluorescent coumarin-benzo[*b*]thiophene 1, 1-dioxide conjugates as mitochondria-targeting antitumor STAT3 inhibitors. *Eur J Med Chem*. 174:236-251.
88. Chen K, Li X, Song G, Zhou T, Long Y, Li Q, Zhong S, Cui Z. Deficiency in the membrane protein Tmbim3a/Grinaa initiates cold-induced ER stress and cell death by activating an intrinsic apoptotic pathway in zebrafish. *J Biol Chem*. 294(30):11445-11457.
89. Xiao Y, Xu D, Song H, Shu F, Wei P, Yang X, Zhong C, Wang X, Müller WE, Zheng Y, Xiao S, Xia Z. Cuprous oxide nanoparticles reduces hypertrophic scarring by inducing fibroblast apoptosis. *INT J NANOMED*. 14:5989-6000.
90. Zhou D, Zhou M, Wang Z, Fu Y, Jia M, Wang X, Liu M, Zhang Y, Sun Y, Lu Y, Tang W, Yi F. PGRN acts as a novel regulator of mitochondrial homeostasis by facilitating mitophagy and mitochondrial biogenesis to prevent podocyte injury in diabetic nephropathy. *Cell Death Dis*. 10(7):524.
91. Li J, Du H, Zhang M, Zhang Z, Teng F, Zhao Y, Zhang W, Yu Y, Feng L, Cui X, Zhang M, Lu T, Guan F, Chen L. Amorphous solid dispersion of Berberine mitigates apoptosis via iPLA2  $\beta$ /Cardiolipin/Opa1 pathway in db/db mice and in Palmitate-treated MIN6  $\beta$ -cells. *Int J Biol Sci*. 15(7):1533-1545.
92. Du L, Lin L, Li Q, Liu K, Huang Y, Wang X, Cao K, Chen X, Cao W, Li F, Shao C, Wang Y, Shi Y. IGF-2 Preprograms Maturing Macrophages to Acquire Oxidative Phosphorylation-Dependent Anti-inflammatory Properties. *Cell Metab*. 29(6):1363-1375.e8.
93. Sun K, Tang S, Hou Y, Xi L, Chen Y, Yin J, Peng M, Zhao M, Cui X, Liu M. Oxidized ATM-mediated glycolysis enhancement in breast cancer-associated fibroblasts contributes to tumor invasion through lactate as metabolic coupling. *EBioMedicine*. 41:370-383.
94. Dai WJ, Qiu J, Sun J, Ma CL, Huang N, Jiang Y, Zeng J, Ren BC, Li WC, Li YH. Downregulation of microRNA-9 reduces inflammatory response and fibroblast proliferation in mice with idiopathic pulmonary fibrosis through the ANO1-mediated TGF- $\beta$ -Smad3 pathway. *J Cell Physiol*. 234(3):2552-2565.
95. Xu B, Wang T, Xiao J, Dong W, Wen HZ, Wang X, Qin Y, Cai N, Zhou Z, Xu J, Wang H. FCPR03, a novel phosphodiesterase 4 inhibitor, alleviates cerebral ischemia/reperfusion injury through activation of the AKT/GSK3 $\beta$ /  $\beta$ -catenin signaling pathway. *Biochem Pharmacol*. 163:234-249.
96. Liu X, Zhao P, Wang X, Wang L, Zhu Y, Gao W. Triptolide Induces Glioma Cell Autophagy and Apoptosis via Upregulating the ROS/JNK and Downregulating the Akt/mTOR Signaling Pathways. *Front Oncol*. 9:387.
97. Wang K, Ma H, Liu H, Ye W, Li Z, Cheng L, Zhang L, Lei Y, Shen L, Zhang F. The Glycoprotein and Nucleocapsid Protein of Hantaviruses Manipulate Autophagy Flux to Restrain Host Innate Immune Responses. *Cell Rep*. 27(7):2075-2091.e5.
98. Yu H, Zhang J, Ji Q, Yu K, Wang P, Song M, Cao Z, Zhang X, Li Y. Melatonin alleviates aluminium chloride-induced immunotoxicity by inhibiting oxidative stress and apoptosis associated with the activation of Nrf2 signaling pathway. *ECOTOX ENVIRON SAFE*. 173:131-141.
99. Shen X, Zhao YF, Xu SQ, Wang L, Cao HM, Cao Y, Zhu Y, Wang Y, Liang ZQ. Cathepsin L induced PC-12 cell apoptosis via activation of B-Myb and regulation of cell cycle proteins. *Acta Pharmacol Sin*. 40(11):1394-1403.
100. Xu WN, Zheng HL, Yang RZ, Liu T, Yu W, Zheng XF, Li B, Jiang SD, Jiang LS. Mitochondrial NDUFA4L2 attenuates the apoptosis of nucleus pulposus cells induced by oxidative stress via the inhibition of mitophagy. *Exp Mol Med*. 51(11):1-16.
101. Zhao W, Xu Z, Cao J, Fu Q, Wu Y, Zhang X, Long Y, Zhang X, Yang Y, Li Y, Mi W. Elamipretide (SS-31) improves mitochondrial dysfunction, synaptic and memory impairment induced by lipopolysaccharide in mice. *J NEUROINFLAMM*. 16(1):230.
102. Sun K, Luo J, Jing X, Guo J, Yao X, Hao X, Ye Y, Liang S, Lin J, Wang G, Guo F. Astaxanthin protects against osteoarthritis via Nrf2: a guardian of cartilage homeostasis. *AGING-US*. 11(22):10513-10531.
103. Fang J, Zhao X, Li S, Xing X, Wang H, Lazarovici P, Zheng W. Protective mechanism of artemisinin on rat bone marrow-derived mesenchymal stem cells against apoptosis induced by hydrogen peroxide via activation of c-Raf-Erk1/2-p90rsk-CREB pathway. *Stem Cell Res Ther*. 10(1):312.
104. Zhang WB, Yang F, Wang Y, Jiao FZ, Zhang HY, Wang LW, Gong ZJ. Inhibition of HDAC6 attenuates LPS-induced inflammation in macrophages by regulating oxidative stress and suppressing the TLR4-MAPK/NF- $\kappa$ B pathways. *Biomed Pharmacother*. 117:109166.
105. Yao W, Lin Z, Wang G, Li S, Chen B, Sui Y, Huang J, Liu Q, Shi P, Lin X, Liu Q, Yao H. Delicaflavone induces apoptosis via mitochondrial pathway accompanying G2/M cycle arrest and inhibition of MAPK signaling cascades in cervical cancer HeLa cells. *Phytomedicine*. 62:152973.
106. Xu WN, Yang RZ, Zheng HL, Yu W, Zheng XF, Li B, Jiang SD, Jiang LS. PGC-1 $\alpha$  acts as a mediator of Sirtuin2 to protect annulus fibrosus from apoptosis induced by oxidative stress through restraining mitophagy. *Int J Biol Macromol*. 136:1007-1017.
107. Xiong S, Sun HJ, Cao L, Zhu M, Liu T, Wu Z, Bian JS. Stimulation of Na<sup>+</sup>/K<sup>+</sup>-ATPase with an Antibody against Its 4th Extracellular Region Attenuates Angiotensin II-Induced H9c2 Cardiomyocyte Hypertrophy via an AMPK/SIRT3/PPAR $\gamma$  Signaling Pathway. *Oxid Med Cell Longev*. 2019:4616034.
108. Yu H, Li JM, Deng K, Zhou W, Wang CX, Wang Q, Li KH, Zhao HY, Huang SW. Tumor acidity activated triphenylphosphonium-based mitochondrial targeting nanocarriers for overcoming drug resistance of cancer therapy. *Theranostics*. 9(23):7033-7050.
109. Wu W, Zhao D, Shah SZA, Zhang X, Lai M, Yang D, Wu X, Guan Z, Li J, Zhao H, Li W, Gao H, Zhou X, Qiao J, Yang L. OPA1 overexpression ameliorates mitochondrial cristae remodeling, mitochondrial dysfunction, and neuronal apoptosis in prion diseases. *Cell Death Dis*. 10(10):710.
110. Yang Q, Sun H, Wang X, Yu X, Zhang J, Guo B, Hexige S. Metabolic changes during malignant transformation in primary cells of oral lichen planus: Succinate accumulation and tumour suppression. *J Cell Mol Med*. 24(2):1179-1188.
111. Dong K, Yang C, Yan Y, Wang P, Sun Y, Wang K, Lu T, Chen Q, Zhang Y, Xing J, Dong Y. Investigation of the intracellular oxidative stress amplification, safety and anti-tumor effect of a kind of novel redox-responsive micelle. *J Mater Chem B*. 6(7):1105-1117.
112. Mengmeng Dong, Jiawen Chen, Jiayuan Zhang, Xiao Liang, Jiebing Yang, Dan Li, Quanshun Li. A Chemoenzymatically Synthesized Cholesterol-G-Poly(amine-Co-Ester)-Mediated p53 Gene Delivery for Achieving Antitumor Efficacy in Prostate Cancer. *INT J NANOMED*. doi: 10.2147/IJN.S191905.
113. Lin Yuan, Xizi Du, Sha Tang, Shuangyan Wu, Leyuan Wang, Yang Xiang, Xiangping Qu, Huijun Liu, Xiaoqun Qin, Chi Liu. ITGB4 Deficiency Induces Senescence of Airway Epithelial Cells Through p53 Activation. *FEBS J*. doi: 10.1111/febs.14749.
114. Jin Wang, Sam Toan, Ruibing Li, Hao Zhou. Melatonin fine-tunes intracellular calcium signals and eliminates myocardial damage through the IP3R/MCU

- pathways in cardiorenal syndrome type 3. *Biochem Pharmacol.* doi: 10.1016/j.bcp.2020.113832.
115. Chong Wang, Mengya Li, Lingling Li, Xiaohui Shen, Yanfang Liu, Shujuan Wang. Stronglyocentrotus nudos Egg Polysaccharide induces autophagy and apoptosis in leukaemia cells by regulating mitochondrial function. *J Cell Mol Med.* doi: 10.1111/jcmm.15995.
116. Xiang Song, Xinzhao Wang, Zhaoyun Liu, Zhiyong Yu. Role of GPX4-Mediated Ferroptosis in the Sensitivity of Triple Negative Breast Cancer Cells to Gefitinib. *Front Oncol.* doi: 10.3389/fonc.2020.597434.
117. Jinqiu Tao, Guangli Sun, Qing Li, Xiaofei Zhi, Zheng Li, Zhongyuan He, Huihui Chen, Aiping Zhou, Jiahui Ye, Guifang Xu, Wenxian Guan, Weijie Zhang. KIF15 promotes the evolution of gastric cancer cells through inhibition of reactive oxygen species-mediated apoptosis. *J Cell Physiol.* doi: 10.1002/jcp.29743.
118. Fang Luo, Sushan Luo, Wenjing Qian, Lin Zhang, Chen Chen, Meimei Xu, Guangling Wang, Zhongfeng Wang, Jian Wang, Wenyuan Wang. Developmental deficits and early signs of neurodegeneration revealed by PD patient derived dopamine neurons. *Stem Cell Res.* doi: 10.1016/j.scr.2020.102027.
119. Xiang Lin, Yongdong Dai, Xiaomei Tong, Wenzhi Xu, Qianmeng Huang, Xiaoying Jin, Chao Li, Feng Zhou, Hanjin Zhou, Xiaona Lin, Dong Huang, Songying Zhang. Excessive oxidative stress in cumulus granulosa cells induced cell senescence contributes to endometriosis-associated infertility. *Redox Biol.* doi: 10.1016/j.redox.2020.101431.
120. Wanyue Huang, Zheng Cao, Qiucheng Yao, Qiang Ji, Jian Zhang, Yanfei Li. Mitochondrial damage are involved in Aflatoxin B 1-induced testicular damage and spermatogenesis disorder in mice. *Sci Total Environ.* doi: 10.1016/j.scitotenv.2019.135077.
121. Wei Li, Dan Li, Hongying Kuang, Xiaoling Feng, Wenxia Ai, Ying Wang, Simao Shi, Jing Chen, Ruifeng Fan. Berberine increases glucose uptake and intracellular ROS levels by promoting Sirtuin 3 ubiquitination. *Biomed Pharmacother.* doi: 10.1016/j.biopha.2019.109563.
122. Dongdong Zhang, Yujiao Liu, Ziyi Luo, Yanling Chen, Anjie Xu, Yuxing Liang, Balu Wu, Xiqin Tong, Xiaoyan Liu, Hui Shen, Li Liu, Yongchang Wei, Haibing Zhou, Yi Liu, Fuling Zhou. The novel thioredoxin reductase inhibitor A-Z2 triggers intrinsic apoptosis and shows efficacy in the treatment of acute myeloid leukemia. *FREE RADICAL BIO MED.* doi: 10.1016/j.freeradbiomed.2019.11.013.
123. Xiaoyan Wang, Zhufeng Chen, Xuemei Fan, Wei Li, Jiaqi Qu, Chuan Dong, Zhixue Wang, Zhenwei Ji, Yang Li. Inhibition of DNMT1L and mitochondrial fission attenuates inflammatory response in fibroblast-like synoviocytes of rheumatoid arthritis. *J Cell Mol Med.* doi: 10.1111/jcmm.14837.
124. Juanjuan Feng, Wenhao Jiang, Yanan Liu, Wanfang Huang, Kewen Hu, Kun Li, Jing Chen, Chengbin Ma, Zhenliang Sun, Xiufeng Pang. Blocking STAT3 by pyrvinium pamoate causes metabolic lethality in KRAS-mutant lung cancer. *Biochem Pharmacol.* doi: 10.1016/j.bcp.2020.113960.
125. Ying Chen, Rui Li, Yuqi Zhu, Sixia Zhong, Jinjun Qian, Dongqing Yang, Artur Jurczynszyn, Meral Beksac, Chunyan Gu, Ye Yang. Dihydroartemisinin Induces Growth Arrest and Overcomes Dexamethasone Resistance in Multiple Myeloma. *Front Oncol.* doi: 10.3389/fonc.2020.00767.
126. Wenjin Guo, Juxiong Liu, Jingxuan Sun, Qian Gong, He Ma, Xingchi Kan, Yu Cao, Jianfa Wang, Shoupeng Fu. Butyrate alleviates oxidative stress by regulating NRF2 nuclear accumulation and H3K9/14 acetylation via GPR109A in bovine mammary epithelial cells and mammary glands. *FREE RADICAL BIO MED.* doi: 10.1016/j.freeradbiomed.2020.01.016.
127. Yunpeng Zhao, Cheng Qiu, Wenhan Wang, Jiangfan Peng, Xiang Cheng, Yangtao Shanguan, Mingyang Xu, Jiayi Li, Ruize Qu, Xiaomin Chen, Suyi Jia, Dan Luo, Long Liu, Peng Li, Fengjin Guo, Krasimir Vasilev, Liang Liu, John Hayball, Shuli Dong, Xin Pan, Yuhua Li, Linlin Guo, Lei Cheng, Weiwei Li. Cortistatin protects against intervertebral disc degeneration through targeting mitochondrial ROS-dependent NLRP3 inflammasome activation. *Theranostics.* doi: 10.7150/thno.45359.
128. Yonglian Zeng, Zhenya Guo, Zhigao Hu, Mingjiang Liu, Yubing Chen, Shilian Chen, Bo Peng, Peng Zhang, Zhan Wu, Hongliang Luo, Fudi Zhong, Keqing Jiang, Yi Lu, Guandou Yuan, Songqing He. FGD1 exhibits oncogenic properties in hepatocellular carcinoma through regulating cell morphology, autophagy and mitochondrial function. *Biomed Pharmacother.* doi: 10.1016/j.biopha.2020.110029.
129. Dejun Xu, Lingbin Liu, Yongju Zhao, Li Yang, Jianyong Cheng, Rongmao Hua, Zelin Zhang, Qingwang Li. Melatonin protects mouse testes from palmitic acid-induced lipotoxicity by attenuating oxidative stress and DNA damage in a SIRT1-dependent manner. *J Pineal Res.* doi: 10.1111/jpi.12690.
130. Min Liu, Xiaoyang Li, Dezhi Huang. Mfn2 Overexpression Attenuates Cardio-Cerebrovascular Ischemia-Reperfusion Injury Through Mitochondrial Fusion and Activation of the AMPK/Sirt3 Signaling. *Front Cell Dev Biol.* doi: 10.3389/fcell.2020.598078.
131. Ranyue Ren, Jiachao Guo, Jia Shi, Yong Tian, Mengwei Li, Hao Kang. PKM2 regulates angiogenesis of VR-EPCs through modulating glycolysis, mitochondrial fission, and fusion. *J Cell Physiol.* doi: 10.1002/jcp.29549.
132. Fang Xu, Mengmeng Guo, Wei Huang, Lili Feng, Jiazhen Zhu, Kangkang Luo, Jian Gao, Bingfeng Zheng, Ling-Dong Kong, Tao Pang, Xudong Wu, Qiang Xu. Annexin A5 regulates hepatic macrophage polarization via directly targeting PKM2 and ameliorates NASH. *Redox Biol.* doi: 10.1016/j.redox.2020.101634.
133. Rui Zhou, Changpei Xiang, Guangzhao Cao, He Xu, Yi Zhang, Hongjun Yang, Jingjing Zhang. Berberine accelerated wound healing by restoring TrxR1/JNK in diabetes. *CLIN SCI.* doi: 10.1042/CS20201145.
134. Wenxing Cui, Xun Wu, Yingwu Shi, Wei Guo, Jianing Luo, Haixiao Liu, Longlong Zheng, Yong Du, Ping Wang, Qiang Wang, Dayun Feng, Shunnan Ge, Yan Qu. 20-HETE synthesis inhibition attenuates traumatic brain injury-induced mitochondrial dysfunction and neuronal apoptosis via the SIRT1/PGC-1  $\alpha$  pathway: A translational study. *CELL PROLIFERAT.* doi: 10.1111/cpr.12964.
135. Xinyu Feng, Shanjie Wang, Xingjun Yang, Jie Lin, Wanrong Man, Yuan Dong, Yan Zhang, Zhijing Zhao, Haichang Wang, Dongdong Sun. Mst1 Knockout Alleviates Mitochondrial Fission and Mitigates Left Ventricular Remodeling in the Development of Diabetic Cardiomyopathy. *Front Cell Dev Biol.* doi: 10.3389/fcell.2020.628842.
136. Doudou Jing, Wei Wu, Xiangyu Deng, Yizhong Peng, Wenbo Yang, Donghua Huang, Zengwu Shao, Dong Zheng. FoxO1a mediated cadmium-induced annulus fibrosus cells apoptosis contributes to intervertebral disc degeneration in smoking. *J Cell Physiol.* doi: 10.1002/jcp.29895.
137. Runze Wang, Yuerong Xu, Wei Zhang, Yexian Fang, Tiquan Yang, Di Zeng, Ting Wei, Jing Liu, Haijia Zhou, Yan Li, Zhan-Peng Huang, Mingming Zhang. Inhibiting miR-22 Alleviates Cardiac Dysfunction by Regulating Sirt1 in Septic Cardiomyopathy. *Front Cell Dev Biol.* doi: 10.3389/fcell.2021.650666.
138. Wanya Li, Botao Liu, Lin Wang, Jilie Liu, Xiuhui Yang, Jia Zheng. Melatonin Attenuates Cardiac Ischemia-Reperfusion Injury through Modulation of IP3R-Mediated Mitochondria-ER Contact. *Oxid Med Cell Longev.* doi: 10.1155/2021/1370862.
139. Jinwen Yuan, Siyan Che, Li Zhang, Xiaomin Li, Junhua Yang, Xiaoming Sun, Zheng Ruan. Assessing the combinatorial cytotoxicity of the exogenous contamination with BDE-209, bisphenol A, and acrylamide via high-content analysis. *Chemosphere.* doi: 10.1016/j.chemosphere.2021.131346.
140. Na Liu, Miao-Miao Lin, Si-Si Huang, Zi-Qi Liu, Jun-Chao Wu, Zhong-Qin Liang, Zheng-Hong Qin, Yan Wang. NADPH and Mito-Apocynin Treatment Protects Against KA-Induced Excitotoxic Injury Through Autophagy Pathway. *Front Cell Dev Biol.* doi: 10.3389/fcell.2021.612554.
141. Changpei Xiang, Fangbo Zhang, Jinhuan Gao, Feifei Guo, Mao Zhang, Rui Zhou, Junying Wei, Ping Wang, Yi Zhang, Jingjing Zhang, Hongjun Yang. Yixin-Shu Capsules Ameliorated Ischemia-Induced Heart Failure by Restoring Trx2 and Inhibiting JNK/p38 Activation. *Oxid Med Cell Longev.* doi: 10.1155/2021/8049079.
142. Xu Wu, Linjing Gong, Liang Xie, Wenyu Gu, Xinyuan Wang, Zilong Liu, Shanqun Li. NLRP3 Deficiency Protects Against Intermittent Hypoxia-Induced Neuroinflammation and Mitochondrial ROS by Promoting the PINK1-Parkin

- Pathway of Mitophagy in a Murine Model of Sleep Apnea. *Front Immunol.* doi: 10.3389/fimmu.2021.628168.
143. Yuhui Liu, Chengjian Shi, Zheng He, Feng Zhu, Min Wang, Ruizhi He, Chunle Zhao, Xiuhui Shi, Min Zhou, Shutao Pan, Yang Gao, Xu Li, Renyi Qin. Inhibition of PI3K/AKT signaling via ROS regulation is involved in Rhein-induced apoptosis and enhancement of oxaliplatin sensitivity in pancreatic cancer cells. *Int J Biol Sci.* doi: 10.7150/ijbs.49514.
144. Lin Zhang, Ruoqi Fu, Dongyu Duan, Ziwei Li, Bin Li, Yue Ming, Li Li, Rui Ni, Jianhong Chen. Cyclovirobuxine D Induces Apoptosis and Mitochondrial Damage in Glioblastoma Cells Through ROS-Mediated Mitochondrial Translocation of Cofilin. *Front Oncol.* doi: 10.3389/fonc.2021.656184.
145. Peng Li, Changlian Xie, Jiankai Zhong, Zhongzhou Guo, Kai Guo, Qiyyun Tu. Melatonin Attenuates ox-LDL-Induced Endothelial Dysfunction by Reducing ER Stress and Inhibiting JNK/Mff Signaling. *Oxid Med Cell Longev.* doi: 10.1155/2021/5589612.
146. Shiyu Song, Shuyin Bao, Chenghong Zhang, Jinwei Zhang, Jiajun Lv, Xiaoyu Li, Maryam Chudhary, Xiang Ren, Li Kong. Stimulation of AMPK Prevents Diabetes-Induced Photoreceptor Cell Degeneration. *Oxid Med Cell Longev.* doi: 10.1155/2021/5587340.
147. Qiang Xie, Jun Zeng, Yongtao Zheng, Tianwen Li, Junwei Ren, Kezhu Chen, Quan Zhang, Rong Xie, Feng Xu, Jianhong Zhu. Mitochondrial Transplantation Attenuates Cerebral Ischemia-Reperfusion Injury: Possible Involvement of Mitochondrial Component Separation. *Oxid Med Cell Longev.* doi: 10.1155/2021/1006636.
148. Fenghe Min, Xin Liu, Yuan Li, Mingyuan Dong, Yidi Qu, Weiwei Liu. Carnosic Acid Suppresses the Development of Oral Squamous Cell Carcinoma via Mitochondrial-Mediated Apoptosis. *Front Oncol.* doi: 10.3389/fonc.2021.760861.
149. Shanshan Dai, Bozhi Ye, Lingfeng Zhong, Yanghao Chen, Guangliang Hong, Guangju Zhao, Lan Su, Zhongqiu Lu. GSDMD Mediates LPS-Induced Septic Myocardial Dysfunction by Regulating ROS-dependent NLRP3 Inflammasome Activation. *Front Cell Dev Biol.* doi: 10.3389/fcell.2021.779432.
150. Rui Huang, Lijun Zhang, Jinmei Jin, Yudong Zhou, Hongwei Zhang, Chao Lv, Dong Lu, Ye Wu, Hong Zhang, Sanhong Liu, Hongzhan Chen, Xin Luan, Weidong Zhang. Bruceine D inhibits HIF-1  $\alpha$ -mediated glucose metabolism in hepatocellular carcinoma by blocking ICAT/  $\beta$ -catenin interaction. *Acta Pharm Sin B.* doi: 10.1016/j.apsb.2021.05.009.
151. Yiyuan Zheng, Chaoyuan Huang, Lina Zhao, Youlan Chen, Fengbin Liu. Regulation of decorin by ursolic acid protects against non-alcoholic steatohepatitis. *Biomed Pharmacother.* doi: 10.1016/j.biopha.2021.112166.
152. Xiaolan Lai, Chuanqian Huang, Xuekun Nie, Qi Chen, Yirong Tang, Xianguo Fu, Ying Lin, Chengjun Nie, Xinyu Xu, Xiukang Wang, Renli Chen, Zichun Chen. Bortezomib Inhibits Multiple Myeloma Cells by Transactivating ATF3 to Trigger miR-135a-5p-Dependent Apoptosis. *Front Oncol.* doi: 10.3389/fonc.2021.720261.
153. Shenglan Zeng, Ze Zhao, Shengnan Zheng, Mengting Wu, Xiaomeng Song, Yiquan Li, Yi Zheng, Bingyu Liu, Lin Chen, Chengjiang Gao, Huiqing Liu. The E3 ubiquitin ligase TRIM31 is involved in cerebral ischemic injury by promoting degradation of TIGAR. *Redox Biol.* doi: 10.1016/j.redox.2021.102058.
154. Shuang Nie, Zhao Shi, Mengyue Shi, Hongzhen Li, Xuettian Qian, Chunyan Peng, Xiwei Ding, Shu Zhang, Ying Lv, Lei Wang, Bo Kong, Xiaoping Zou, Shanshan Shen. PPAR $\gamma$ /SOD2 Protects Against Mitochondrial ROS-Dependent Apoptosis via Inhibiting ATG4D-Mediated Mitophagy to Promote Pancreatic Cancer Proliferation. *Front Cell Dev Biol.* doi: 10.3389/fcell.2021.745554.
155. Wen-Jun Liu, Peng-Yu Pan, Ye Sun, Jian-Bo Wang, Huan Zhou, Xin Xie, Zhi-Yuan Duan, Han-Yu Dong, Wen-Na Chen, Li-de Zhang, Chun Wang. Deferoxamine Counteracts Cisplatin Resistance in A549 Lung Adenocarcinoma Cells by Increasing Vulnerability to Glutamine Deprivation-Induced Cell Death. *Front Oncol.* doi: 10.3389/fonc.2021.794735.
156. Congcong Wang, Lei Wang, Chengmeng Huang, Yungang Liu, Jian Liu, Hongxuan Kuang, Qihua Pang, Hongyu Han, Ruifang Fan. Involvement of NLRP3/Caspase-1/GSDMD-Dependent pyroptosis in BPA-Induced apoptosis of human neuroblastoma cells. *Biochem Pharmacol.* doi: 10.1016/j.bcp.2022.115042.
157. Liang Dong, Yun-Jun Xu, Cong Sui, Yang Zhao, Li-Bo Mao, Denis Gebauer, Rose Rosenberg, Jonathan Avaro, Ya-Dong Wu, Huai-Ling Gao, Zhao Pan, Hui-Qin Wen, Xu Yan, Fei Li, Yang Lu, Helmut Cölfen, Shu-Hong Yu. Highly hydrated paramagnetic amorphous calcium carbonate nanoclusters as an MRI contrast agent. *Nat Commun.* doi: 10.1038/s41467-022-32615-3.
158. Kuan Liu, Zhiqi Liu, Zhuofan Liu, Zhuo Ma, Yu Deng, Wei Liu, Bin Xu. Manganese induces S-nitrosylation of PINK1 leading to nerve cell damage by repressing PINK1/Parkin-mediated mitophagy. *Sci Total Environ.* doi: 10.1016/j.scitotenv.2022.155358.
159. Yuqi Zhang, Enze Yao, Yijing Liu, Yining Zhang, Mengyang Ding, Jingyu Liu, Xiaoming Chen, Sairong Fan. FUT2 Facilitates Autophagy and Suppresses Apoptosis via p53 and JNK Signaling in Lung Adenocarcinoma Cells. *Cells.* doi: 10.3390/cells11244031.
160. Ruzhou Zhao, Yixin Xu, Xiaobo Wang, Xiang Zhou, Yanqi Liu, Shuai Jiang, Lin Zhang, Zhibin Yu. Withaferin A Enhances Mitochondrial Biogenesis and BNIP3-Mediated Mitophagy to Promote Rapid Adaptation to Extreme Hypoxia. *Cells.* doi: 10.3390/cells12010085.
161. Dandan Liu, Jiande Li, Xiaoming Rong, Jie Li, Ying Peng, Qingyu Shen. Cdk5 Promotes Mitochondrial Fission via Drp1 Phosphorylation at S616 in Chronic Ethanol Exposure-Induced Cognitive Impairment. *Mol Neurobiol.* doi: 10.1007/s12035-022-03008-w.
162. Fei Jia, Yu Liu, Xinyu Dou, Chuanchao Du, Tianli Mao, Xiaoguang Liu. Liensinine Inhibits Osteosarcoma Growth by ROS-Mediated Suppression of the JAK2/STAT3 Signaling Pathway. *Oxid Med Cell Longev.* doi: 10.1155/2022/8245614.
163. Bonan Xiao, Yilong Cui, Bo Li, Jian Zhang, Xuliang Zhang, Miao Song, Yanfei Li. ROS antagonizes the protection of Parkin-mediated mitophagy against aluminum-induced liver inflammatory injury in mice. *Food Chem Toxicol.* doi: 10.1016/j.fct.2022.113126.
164. Qi Wang, Fubo Jia, Chen Guo, Yuping Wang, Xuliang Zhang, Yilong Cui, Miao Song, Zheng Cao, Yanfei Li. PINK1/Parkin-mediated mitophagy as a protective mechanism against AFB1-induced liver injury in mice. *Food Chem Toxicol.* doi: 10.1016/j.fct.2022.113043.
165. Yan-Ran Sheng, Wen-Ting Hu, Hui-Hui Shen, Chun-Yan Wei, Yu-Kai Liu, Xiao-Qian Ma, Ming-Qing Li, Xiao-Yong Zhu. An imbalance of the IL-33/ST2-AXL-efferocytosis axis induces pregnancy loss through metabolic reprogramming of decidual macrophages. *Cell Mol Life Sci.* doi: 10.1007/s00018-022-04197-2.
166. Meng Duan, Hainan Chen, Linjie Yin, Xiao Zhu, Petr Novák, Yuncheng Lv, Guojun Zhao, Kai Yin. Mitochondrial apolipoprotein A-I binding protein alleviates atherosclerosis by regulating mitophagy and macrophage polarization. *Cell Commun Signal.* doi: 10.1186/s12964-022-00858-8.
167. Liling Wang, Huan Wang, Jing Luo, Ting Xie, Gil Mor, Aihua Liao. Decorin promotes decidual M1-like macrophage polarization via mitochondrial dysfunction resulting in recurrent pregnancy loss. *Theranostics.* doi: 10.7150/thno.78467.
168. Zhitao Jiang, Xiang Yang, Yi Han, Jie Li, Chen Hu, Chundi Liu, Wei Xiao. Sarmenosin promotes USP17 and regulates Nrf2-mediated mitophagy and cellular oxidative stress to alleviate APAP-induced acute liver failure. *Phytomedicine.* doi: 10.1016/j.phymed.2022.154337.

注：更多使用本产品的文献请参考产品网页。

Version 2024.03.20