碧云天生物技术/Beyotime Biotechnology  
订货热线: 400-1683301或800-8283301  
订货e-mail: order@beyotime.com  
技术咨询: info@beyotime.com  
网址: http://www.beyotime.com

## AdPlus-mCherry-GFP-LC3B

产品编号	产品名称	包装
C3012-1ml	AdPlus-mCherry-GFP-LC3B	1ml
C3012-10ml	AdPlus-mCherry-GFP-LC3B	10ml

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Table 1. Summary of Recommended Biosafety Levels for Infectious Agents

BSL	Agents	Practices	Primary Barriers and Safety Equipment	Facilities (Secondary Barriers)
1	Not known to consistently cause diseases in healthy adults	Standard microbiological practices	<ul style="list-style-type: none"> <li>■ No primary barriers required.</li> <li>■ PPE: laboratory coats and gloves; eye, face protection, as needed</li> </ul>	Laboratory bench and sink required
2	<ul style="list-style-type: none"> <li>■ Agents associated with human disease</li> <li>■ Routes of transmission include percutaneous injury, ingestion, mucous membrane exposure</li> </ul>	BSL-1 practice plus: <ul style="list-style-type: none"> <li>■ Limited access</li> <li>■ Biohazard warning signs</li> <li>■ “Sharps” precautions</li> <li>■ Biosafety manual defining any needed waste decontamination or medical surveillance policies</li> </ul>	Primary barriers: <ul style="list-style-type: none"> <li>■ BSCs or other physical containment devices used for all manipulations of agents that cause splashes or aerosols of infectious materials</li> <li>■ PPE: Laboratory coats, gloves, face and eye protection, as needed</li> </ul>	BSL-1 plus: <ul style="list-style-type: none"> <li>■ Autoclave available</li> </ul>
3	Indigenous or exotic agents that may cause serious or potentially lethal disease through the inhalation route of exposure	BSL-2 practice plus: <ul style="list-style-type: none"> <li>■ Controlled access</li> <li>■ Decontamination of all waste</li> <li>■ Decontamination of laboratory clothing before laundering</li> </ul>	Primary barriers: <ul style="list-style-type: none"> <li>■ BSCs or other physical containment devices used for all open manipulations of agents</li> <li>■ PPE: Protective laboratory clothing, gloves, face, eye and respiratory protection, as needed</li> </ul>	BSL-2 plus: <ul style="list-style-type: none"> <li>■ Physical separation from access corridors</li> <li>■ Self-closing, double-door access</li> <li>■ Exhausted air not recirculated</li> <li>■ Negative airflow into laboratory</li> <li>■ Entry through airlock or anteroom</li> <li>■ Hand washing sink near laboratory exit</li> </ul>

<p><b>4</b> ■ Dangerous/exotic agents which post high individual risk of aerosol-transmitted laboratory infections that are frequently fatal, for which there are no vaccines or treatments</p> <ul style="list-style-type: none"> <li>■ Agents with a close or identical antigenic relationship to an agent requiring BSL-4 until data are available to redesignate the level</li> <li>■ Related agents with unknown risk of transmission</li> </ul>	<p>BSL-3 practices plus:</p> <ul style="list-style-type: none"> <li>■ Clothing change before entering</li> <li>■ Shower on exit</li> <li>■ All material decontaminated on exit from facility</li> </ul>	<p>Primary barriers:</p> <ul style="list-style-type: none"> <li>■ All procedures conducted in Class III BSCs or Class I or II BSCs in combination with full-body, air-supplied, positive pressure suit</li> </ul>	<p>BSL-3 plus:</p> <ul style="list-style-type: none"> <li>■ Separate building or isolated zone</li> <li>■ Dedicated supply and exhaust, vacuum, and decontamination systems</li> <li>■ Other requirements outlined in the text</li> </ul>
---	---	---	---

BSL, biosafety level; PPE, personal protective equipment.

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产品编号	产品名称	包装
C3006-1ml	Ad-GFP-LC3B	1ml
C3006-10ml	Ad-GFP-LC3B	10ml
C3011-1ml	Ad-mCherry-GFP-LC3B	1ml
C3011-10ml	Ad-mCherry-GFP-LC3B	10ml
C3012-1ml	AdPlus-mCherry-GFP-LC3B	1ml
C3012-10ml	AdPlus-mCherry-GFP-LC3B	10ml

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1. Yiyu Cai, Xinzhu Li, Zihao Pan, Yue Zhu, Ji Tuo, Qinghai Meng, Guoliang Dai, Guangming Yang, Yang Pan. Anthocyanin ameliorates hypoxia and ischemia induced inflammation and apoptosis by increasing autophagic flux in SH-SY5Y cells Eur J Pharmacol. 2020 Sep 15;883:173360.;doi: 10.1016/j.ejphar.2020.173360.

Version 2021.09.01