



碧云天生物技术/Beyotime Biotechnology
订货热线: 400-1683301 或 800-8283301
订货 e-mail: order@beyotime.com
技术咨询: info@beyotime.com
网址: <http://www.beyotime.com>

RIPA裂解液(强)

产品编号	产品名称	包装
P0013B	RIPA裂解液(强)	100ml

产品简介:

- 碧云天生产的RIPA裂解液(RIPA Lysis Buffer)是一种传统的细胞组织快速裂解液。RIPA裂解液裂解得到的蛋白样品可以用于常规的PAGE、Western、免疫沉淀(immunoprecipitation, IP)、免疫共沉淀(co-IP)和ELISA等。
- 本产品可以用于动物、植物的细胞或组织样品,也可以用于真菌或细菌样品。
- RIPA的本意是Radio Immunoprecipitation Assay。RIPA裂解液的配方有很多种,根据其裂解液的强度大致可以分为强、中、弱三类。关于不同的RIPA裂解液以及碧云天生产的其它裂解液的主要特点和差异,以及如何选择裂解液可参考我们的相关网页:<http://www.beyotime.com/support/lysis-buffer.htm>。
- RIPA裂解液(强)的主要成分为50mM Tris (pH 7.4), 150mM NaCl, 1% Triton X-100, 1% sodium deoxycholate, 0.1% SDS, 以及sodium orthovanadate, sodium fluoride, EDTA, leupeptin等多种抑制剂。可以有效抑制蛋白降解。
- 用RIPA裂解液裂解得到的蛋白样品,可以用碧云天生产的BCA蛋白浓度测定试剂盒(P0009/P0010/P0010S/P0011/P0012/P0012S)测定蛋白浓度。由于含有较高浓度的去垢剂,不能用Bradford法测定由本裂解液裂解得到样品的蛋白浓度。

包装清单:

产品编号	产品名称	包装
P0013B	RIPA裂解液(强)	100ml
—	说明书	1份

保存条件:

-20°C保存,一年有效。

注意事项:

- 为取得最佳的使用效果,尽量避免过多的反复冻融。可以适当分装后使用。
- 需自备PMSF。PMSF(ST506)可以向碧云天订购。也可以选购总体效果更佳的碧云天生产的P1045/P1046 蛋白酶磷酸酶抑制剂混合物(通用型, 50X), 或者根据具体用途选择P1048/P1049 蛋白酶磷酸酶抑制剂混合物(通用型, 质谱兼容, 50X)、P1050/P1051蛋白酶磷酸酶抑制剂混合物(哺乳动物样品抽提用, 50X)、P1055/P1056 蛋白酶磷酸酶抑制剂混合物(植物样品抽提用, 50X)、P1060/P1061 蛋白酶磷酸酶抑制剂混合物(真菌或酵母抽提用, 50X)、P1065/P1066 蛋白酶磷酸酶抑制剂混合物(细菌抽提用, 50X)。如果无需检测磷酸化蛋白, 也可以选不含磷酸酶抑制剂的蛋白酶抑制剂化合物。
- 裂解样品的所有步骤都需在冰上或4°C进行。
- 关于裂解液的选择, 一方面可以参考碧云天的相关网页: <http://www.beyotime.com/support/lysis-buffer.htm>选择合适的裂解液; 另一方面也需要通过一些预实验来摸索最佳的适合您实验条件的裂解液。
- 本产品仅限于专业人员的科学研究用, 不得用于临床诊断或治疗, 不得用于食品或药品, 不得存放于普通住宅内。
- 为了您的安全和健康, 请穿实验服并戴一次性手套操作。

使用说明:

1. 对于培养细胞样品:

- 融解RIPA裂解液, 混匀。取适当量的裂解液, 在使用前数分钟内加入PMSF, 使PMSF的最终浓度为1mM, 或者根据实验需要加入适当的上述蛋白酶磷酸酶抑制剂混合物。
- 对于贴壁细胞: 去除培养液, 用PBS、生理盐水或无血清培养液洗一遍(如果血清中的蛋白没有干扰, 可以不洗)。按照6孔板每孔加入150-250微升裂解液的比例加入裂解液。用枪吹打数下, 使裂解液和细胞充分接触。通常裂解液接触动物细胞1-2秒后, 细胞就会被裂解。植物细胞宜在冰上裂解2-10min。
对于悬浮细胞: 离心收集细胞, 轻轻vortex或者弹击管底以把细胞尽量分散开。按照6孔板每孔细胞加入150-250微升裂解液的比例加入裂解液。轻弹管底以充分裂解细胞。充分裂解后应没有明显的细胞沉淀。如果细胞量较多, 必需分装成50-100万细胞/管, 然后再裂解。
对于细菌或酵母: 对于1ml菌液或酵母液, 离心去上清, 如果有必要可以使用PBS洗涤一次, 充分去除液体后, 轻轻vortex或者弹击管底以把细菌或酵母尽量弹散。加入100-200微升裂解液, 轻轻vortex或者弹击管底以混匀, 冰上裂解2-10min。如果希望获得更好的裂解效果, 细菌和酵母可以分别使用溶菌酶和破壁酶(Lyticase)消化, 然后再使用本裂解液进行裂解。

裂解液用量说明：通常6孔板每孔细胞或者1ml的菌液或酵母液中的细菌和酵母量加入150微升裂解液已经足够，但如果细胞密度非常高可以适当加大裂解液的用量到200微升或250微升。每100万动物细胞用100微升本产品裂解后获得的上清，其蛋白浓度约为2-4mg/ml，不同细胞有所不同。

c. 充分裂解后，10000-14000g离心3-5分钟，取上清，即可进行后续的PAGE、Western和免疫沉淀等操作。

2. 对于组织样品：

a. 把组织剪切成细小的碎片。

b. 融解RIPA裂解液，混匀。取适当量的裂解液，在使用前数分钟内加入PMSF，使PMSF的最终浓度为1mM，或者根据实验需要加入适当的上述蛋白酶抑制剂混合物。

c. 按照每20毫克组织加入150-250微升裂解液的比例加入裂解液。(如果裂解不充分可以适当添加更多的裂解液，如果需要高浓度的蛋白样品，可以适当减少裂解液的用量。)

d. 用玻璃匀浆器匀浆，或使用碧云天生产的E6600 TissueMaster™手持式组织研磨仪研磨，直至充分裂解。也可以把组织样品冷冻后液氮研磨，研磨充分后加入裂解液进行裂解。

e. 充分裂解后，10000-14000g离心3-5分钟，取上清，即可进行后续的PAGE、Western和免疫沉淀等操作。每20mg冻存的小鼠肝脏组织用200微升本裂解液裂解后获得的上清，其蛋白浓度约为15-25mg/ml，不同状态的不同组织有所不同。

f. 如果组织样品本身非常细小，可以适当剪切后直接加入裂解液裂解，通过强烈vortex使样品裂解充分。然后同样离心取上清，用于后续实验。直接裂解的优点是比较方便，不必使用匀浆器或研磨设备，缺点是不如匀浆或研磨那样裂解得比较充分。

注：RIPA裂解液的裂解产物中经常会出现一小团透明胶状物，属正常现象。该透明胶状物为含有基因组DNA等的复合物。在不检测和基因组DNA结合特别紧密的蛋白的情况下，可以直接离心取上清用于后续实验；如果需要检测和基因组结合特别紧密的蛋白，则可以通过超声处理打碎打散该透明胶状物，随后离心取上清用于后续实验。如果检测一些常见的转录因子，例如NF-kappaB、p53等时，通常不必进行超声处理，就可以检测到这些转录因子。

附录：碧云天生产的各种裂解液主要特点、差异和选择

首先请参考下表，了解各种裂解液的主要特点和差异。

产品编号	P0013	P0013B	P0013C	P0013D	P0013F	P0013G	P0013J	P0013K
产品名称	Western及IP细胞裂解液	RIPA裂解液(强)	RIPA裂解液(中)	RIPA裂解液(弱)	NP-40裂解液	SDS裂解液	Western及IP细胞裂解液(无抑制剂)	RIPA裂解液(强, 无抑制剂)
有效裂解成分	1% Triton X-100	1% Triton X-100, 1% deoxycholate, 0.1% SDS	1% NP-40, 0.5% deoxycholate, 0.1% SDS	1% NP-40, 0.25% deoxycholate	1% NP-40	1% SDS	1% Triton X-100	1% Triton X-100, 1% deoxycholate, 0.1% SDS
裂解强度	温和	强	中	温和	温和	强	温和	强
对膜蛋白的提取	一般	很好	较好	一般	一般	很好	一般	很好
对胞浆蛋白的提取	很好	很好	很好	很好	很好	很好	很好	很好
对核蛋白的提取	较好	很好	较好	较好	较好	很好	较好	很好
胞浆磷酸化蛋白提取	很好	很好	很好	很好	很好	很好	很好	很好
细胞核转录因子提取	很好	很好	很好	很好	很好	很好	很好	很好
含蛋白酶抑制剂	是	是	是	是	是	是	否	否
含磷酸酯酶抑制剂	是	是	是	是	是	是	否	否
不同物种样品兼容性	高	高	高	高	高	高	高	高
主要用途	WB, IP, co-IP	WB, IP	WB, IP	WB, IP, co-IP	WB, IP, co-IP	WB, ChIP	WB, IP, co-IP	WB, IP

➤ 用于普通的Western、IP或co-IP，我们推荐使用Western及IP细胞裂解液(P0013)，该裂解液已被国内各大研究机构广泛使用，发表大量SCI论文，用户普遍反映很好。裂解细胞或组织后，没有非常粘滞的透明状DNA团块形成，不必采用超声处理等就可以非常理想地用于后续操作。另外该裂解液裂解的产物也适合用于磷酸化蛋白的Western检测。

➤ 对于某些特殊蛋白的IP，如果发现Western及IP细胞裂解液(P0013)效果不是非常理想，可以尝试用RIPA裂解液(强、中或弱)或NP-40裂解液。如果发现IP的时候背景很高，即非特异的蛋白也被IP下来，则需要选用裂解强度较高的裂解液，例如RIPA裂解液(强或中)。如果发现目的蛋白无法被IP下来，则说明裂解液的强度过强，可以使用较为温和的裂解液例如RIPA裂解液(弱)或NP-40裂解液。

➤ 对于某些难溶解蛋白的Western，如果发现Western及IP细胞裂解液(P0013)效果不是非常理想，可以尝试使用裂解强度更高的裂解液例如RIPA裂解液(强、中)或SDS裂解液。使用RIPA裂解液(强)的用户也非常多，发表了大量SCI论文。

➤ 用于特定用途需要自行添加特定抑制剂或不需要添加抑制剂时，可以考虑选购P0013J或P0013K。P0013J在很多时候可以兼容酶活性和生物小分子的检测，对于特定的酶或生物小分子的检测是否兼容需要自行测试，碧云天不提供具体的应用信息。P0013J的裂解能力比P0013K弱一些，但用于酶活性和生物小分子时，P0013J的兼容性通常会更好一些。

使用本产品的文献:

- Deng XQ, Chen LL, Li NX. The expression of SIRT1 in nonalcoholic fatty liver disease induced by high-fat diet in rats. *Liver Int* 2007 Jun;27(5):708-15.
- Shen Y, Yang X, Dong N, Xie X, Bai X, Shi Y. Generation and selection of immunized Fab phage display library against human B cell lymphoma. *Cell Res* 2007 Jul;17(7):650-60.
- Qin ZX, Zhu HY, Hu YH. Effects of lysophosphatidylcholine on beta-amyloid-induced neuronal apoptosis. *Acta Pharmacol Sin* 2009 Apr;30(4):388-95.
- Chen Y, Zhang H, Gou X, Horikawa Y, Xing J, Chen Z. Upregulation of HAB18G/CD147 in activated human umbilical vein endothelial cells enhances the angiogenesis. *Cancer Lett* 2009 Jun 8;278(1):113-21.
- Liu RT, Zou LB, Lü QJ. Liquiritigenin inhibits Abeta(25-35)-induced neurotoxicity and secretion of Abeta(1-40) in rat hippocampal neurons. *Acta Pharmacol Sin* 2009 Jul;30(7):899-906.
- Xiao X, Liu A, Wen H, Tian Y, Ni J, Liu G. Expression and localization of transcription factor Ets-1 in the rat ovary during the estrous cycle and pregnancy. *Fertil Steril* 2009;91(5 Suppl):1998-2005.
- Chen W, Hou J, Yin Y, Jang J, Zheng Z, Fan H, Zou G. alpha-Bisabolol induces dose- and time-dependent apoptosis in HepG2 cells via a Fas- and mitochondrial-related pathway, involves p53 and NFkappaB. *Biochem Pharmacol* 2010;80(2):247-54.
- Li XC, Tong GX, Zhang Y, Liu SX, Jin QH, Chen HH, Chen P. Neferine inhibits angiotensin II-stimulated proliferation in vascular smooth muscle cells through heme oxygenase-1. *Acta Pharmacol Sin* 2010;31(6):679-86.
- Wu W, Yan C, Gan T, Chen Z, Lu X, Duerksen-Hughes PJ, Zhu X, Yang J. Nuclear proteome analysis of cisplatin-treated HeLa cells. *Mutat Res-Rev Mutat* 2010;691(1-2):1-8.
- J. Dang, R.Jia, Y.Tu, S.Xiao, G.Ding. Erythropoietin prevents reactive oxygen species generation and renal tubular cell apoptosis at high glucose level. *Biomed Pharmacother* 2010 Dec;64(10):681-5.
- Ma Y, Niu R, Sun Z, Wang J, Luo G, Zhang J, Wang J. Inflammatory responses induced by fluoride and arsenic at toxic concentration in rabbit aorta. *Arch Toxicol* 2012 Jun;86(6):849-56.
- Yu X, Tang J, Wang Q, Ye W, Tao K, Duan S, Lu C, Yang X, Dong S, Zheng X, Wang Y. The RxLR effector Avh241 from *Phytophthora sojae* requires plasma membrane localization to induce plant cell death. *New Phytol* 2012 Oct;196(1):247-60.
- Wang Y, Liu Q, Liu Z, Li B, Sun Z, Zhou H, Zhang X, Gong Y, Shao C. Berberine, a genotoxic alkaloid, induces ATM-Chk1 mediated G2 arrest in prostate cancer cells. *Mutat Res-Rev Mutat* 2012 Jun 1;734(1-2):20-9.
- Li R, Yu C, Gao R, Liu X, Lu J, Zhao L, Chen X, Ding Y, Wang Y, He J. Effects of DEHP on endometrial receptivity and embryo implantation in pregnant mice. *J Hazard Mater* 2012 Nov 30;241-242:231-40.
- Gao L, Song JR, Zhang JW, Zhao X, Zhao QD, Sun K, Deng WJ, Li R, Lv G, Cheng HY, Wei LX. Chloroquine promotes the anticancer effect of TACE in a rabbit VX2 liver tumor model. *Int J Biol Sci* 2013;9(4):322-30.
- Pang X, Wei Y, Zhang Y, Zhang M, Lu Y, Shen P. Peroxisome proliferator-activated receptor- γ activation inhibits hepatocellular carcinoma cell invasion by upregulating plasminogen activator inhibitor-1. *Cancer Sci* 2013 Jun;104(6):672-80.
- Xu P, Zhang XX, Miao C, Fu Z, Li Z, Zhang G, Zheng M, Liu Y, Yang L, Wang T. Promotion of melanoma cell invasion and tumor metastasis by microcystin-LR via phosphatidylinositol 3-kinase/AKT pathway. *Environ Sci Technol* 2013 Aug 6;47(15):8801-8.
- Mo QQ, Chen PB, Jin X, Chen Q, Tang L, Wang BB, Li KZ, Wu P, Fang Y, Wang SX, Zhou JF, Ma D, Chen G. Inhibition of Hecl expression enhances the sensitivity of human ovarian cancer cells to paclitaxel. *Acta Pharmacol Sin* 2013 Apr;34(4):541-8.
- Liu E, Kuang Y, He W, Xing X, Gu J. Casticin Induces Human Glioma Cell Death through Apoptosis and Mitotic Arrest. *CELL PHYSIOL BIOCHEM* 2013 Jun 3;31(6):805-814.
- Xu WP, Yi M, Li QQ, Zhou WP, Cong WM, Yang Y, Ning BF, Yin C, Huang ZW, Wang J, Qian H, Jiang CF, Chen YX, Xia CY, Wang HY, Zhang X, Xie WF. Perturbation of miR-370-LIN28A-NF- κ B regulatory circuit contributes to the development of hepatocellular carcinoma. *Hepatology* 2013 Dec;58(6):1977-91.
- Yin C, Wang PQ, Xu WP, Yang Y, Zhang Q, Ning BF, Zhang PP, Zhou WP, Xie WF, Chen WS, Zhang X. Hepatocyte nuclear factor-4a reverses malignancy of hepatocellular carcinoma through regulating miR-134 in the DLK1-DIO3 region. *Hepatology* 2013 Dec;58(6):1964-76.
- Zhang X, Chen G, Lu Y, Liu J, Fang M, Luo J, Cao Q, Wang X. Association of Mitochondrial Letm1 with Epileptic Seizures. *Cereb Cortex* 2014 Oct;24(10):2533-40.
- Sun H, Wu Y, Fu D, Liu Y, Huang C. SIRT6 regulates osteogenic differentiation of rat bone marrow mesenchymal stem cells partially via suppressing the nuclear factor- κ B signaling pathway. *Stem Cells* 2014 Jul;32(7):1943-55.
- Cao M, Hou D, Liang H, Gong F, Wang Y, Yan X, Jiang X, Wang C, Zhang J, Zen K, Zhang CY, Chen X. miR-150 promotes the proliferation and migration of lung cancer cells by targeting SRC kinase signalling inhibitor 1. *Eur J Cancer* 2014 Mar;50(5):1013-24.
- Wang G, Liu X, Meng L, Liu S, Wang L, Li J, Cui C, Meng J, Hu S, Wei Y. Up-regulated lipocalin-2 in pulmonary hypertension involving in pulmonary artery SMC resistance to apoptosis. *Int J Biol Sci* 2014 Jul 17;10(7):798-806.
- Deng Z, Yan F, Jin Q, Li F, Wu J, Liu X, Zheng H. Reversal of multidrug resistance phenotype in human breast cancer cells using doxorubicin-liposome-microbubble complexes assisted by ultrasound. *J Control Release* 2014 Jan 28;174:109-16.
- Guo S, Xu X, Tang Y, Zhang C, Li J, Ouyang Y, Ju J, Bie P, Wang H. miR-15a inhibits cell proliferation and epithelial to mesenchymal transition in pancreatic ductal adenocarcinoma by down-regulating Bmi-1 expression. *Cancer Lett* 2014 Mar 1;344(1):40-6.
- Cong ZX, Wang HD, Zhou Y, Wang JW, Pan H, Zhang DD, Zhang L, Zhu L. Temozolomide and irradiation combined treatment-induced Nrf2 activation increases chemoradiation sensitivity in human glioblastoma cells. *Neuro-Oncology* 2014 Jan;116(1):41-8.
- Wang S, Li Y, Fan J, Wang Z, Zeng X, Sun Y, Song P, Ju D. The role of autophagy in the neurotoxicity of cationic PAMAM dendrimers. *Biomaterials* 2014 Aug;35(26):7588-97.
- Zhang Y, Lv X, Bai Y, Zhu X, Wu X, Chao J, Duan M, Buch S, Chen L, Yao H. Involvement of sigma-1 receptor in astrocyte activation induced by methamphetamine via up-regulation of its own expression. *J Neuroinflamm* 2015 Feb 17;12:29.
- Xie Q, Wang Z, Huang Y, Bi X, Zhou H, Lin M, Yu Z, Wang Y, Ni N, Sun J, Wu S, You Z, Guo C, Sun H, Wang Y, Gu P, Fan X. Characterization of human ethmoid sinus mucosa derived mesenchymal stem cells (hESMSCs) and the application of hESMSCs cell sheets in bone regeneration. *Biomaterials* 2015 Oct;66:67-82.
- Liu GY, Jiang XX, Zhu X, He WY, Kuang YL, Ren K, Lin Y, Gou X. ROS activates JNK-mediated autophagy to counteract apoptosis in mouse mesenchymal stem cells in vitro. *Acta Pharmacol Sin* 2015 Dec;36(12):1473-9.
- Xie Q, Wang Z, Zhou H, Yu Z, Huang Y, Sun H, Bi X, Wang Y, Shi W, Gu P, Fan X. The role of miR-135-modified adipose-derived mesenchymal stem cells in bone regeneration. *Biomaterials* 2016 Jan;75:279-94.
- Wang Q, Zhang X, Shen E, Gao J, Cao F, Wang X, Li Y, Tian T, Wang J, Chen Z, Wang J, Shen L. The anti-HER3 antibody in combination with trastuzumab exerts synergistic antitumor activity in HER2-positive gastric cancer. *Cancer Lett* 2016 Sep 28;380(1):20-30.
- Cheng X, He S, Yuan J, Miao S, Gao H, Zhang J, Li Y, Peng W, Wu P. Lipoxin A4 attenuates LPS-induced mouse acute lung injury via Nrf2-mediated E-cadherin expression in airway epithelial cells. *Free Radical Bio Med* 2016 Apr;93:52-66.
- Wang WT, Ye H, Wei PP, Han BW, He B, Chen ZH, Chen YQ. LncRNAs H19 and

- HULC, activated by oxidative stress, promote cell migration and invasion in cholangiocarcinoma through a ceRNA manner. *J Hematol Oncol* 2016 Nov 3;9(1):117.
37. Min J, Weitian Z, Peng C, Yan P, Bo Z, Yan W, Yun B, Xukai W. Correlation between insulin-induced estrogen receptor methylation and atherosclerosis. *Cardiovasc Diabetol* 2016 Nov 10;15(1):156.
 38. Qu Z, Wu J, Wu J, Luo D, Jiang C, Ding Y. Exosomes derived from HCC cells induce sorafenib resistance in hepatocellular carcinoma both in vivo and in vitro. *J Exp Clin Canc Res* 2016 Sep 30;35(1):159.
 39. Chen Y, Zhao R, Zhao Q, Shao Y, Zhang S. Knockdown of HPIP Inhibits the Proliferation and Invasion of Head-and-Neck Squamous Cell Carcinoma Cells by Regulating PI3K/Akt Signaling Pathway. *Oncol Res* 2016;24(3):153-60.
 40. Liu L, Li Y, Wang Y, Zhao P, Wei S, Li Z, Chang H, He H. Biochemical characterization and functional analysis of the POU transcription factor POU-M2 of *Bombyx mori*. *Int J Biol Macromol* 2016 May;86:701-8.
 41. Liu W, Zhang Y, Xia P, Li S, Feng X, Gao Y, Wang K, Song Y, Duan Z, Yang S, Shao Z, Yang C. MicroRNA-7 regulates IL-1 β -induced extracellular matrix degeneration by targeting GDF5 in human nucleus pulposus cells. *Biomed Pharmacother* 2016 Oct;83:1414-1421.
 42. Lu Y, Xu S, Chen H, He M, Deng Y, Cao Z, Pi H, Chen C, Li M, Ma Q, Gao P, Ji Y, Zhang L, Yu Z, Zhou Z. CdSe/ZnS quantum dots induce hepatocyte pyroptosis and liver inflammation via NLRP3 inflammasome activation. *Biomaterials* 2016 Jun;90:27-39.
 43. Fan TF, Wu TF, Bu LL, Ma SR, Li YC, Mao L, Sun ZJ, Zhang WF. Dihydromyricetin promotes autophagy and apoptosis through ROS-STAT3 signaling in head and neck squamous cell carcinoma. *ONCOTARGET* 2016 Sep 13;7(37):59691-59703.
 44. Shen ZL, Wang B, Jiang KW, Ye CX, Cheng C, Yan YC, Zhang JZ, Yang Y, Gao ZD, Ye YJ, Wang S. Downregulation of miR-199b is associated with distant metastasis in colorectal cancer via activation of SIRT1 and inhibition of CREB/KISS1 signaling. *ONCOTARGET* 2016 Jun 7;7(23):35092-105.
 45. Wang Y, Leng H, Chen H, Wang L, Jiang N, Huo X, Yu B. Knockdown of UBE2T Inhibits Osteosarcoma Cell Proliferation, Migration, and Invasion by Suppressing the PI3K/Akt Signaling Pathway. *Oncol Res* 2016;24(5):361-369.
 46. Xie S, Jiang H, Zhai XW, Wei F, Wang SD, Ding J, Chen Y. Antitumor action of CDK inhibitor LS-007 as a single agent and in combination with ABT-199 against human acute leukemia cells. *Acta Pharmacol Sin* 2016 Nov;37(11):1481-1489.
 47. Zhao Z, Zhao J, Xue J, Zhao X, Liu P. Autophagy inhibition promotes epithelial-mesenchymal transition through ROS/HO-1 pathway in ovarian cancer cells. *Am J Cancer Res* 2016 Oct 1;6(10):2162-2177.
 48. Huang W, Li J, Guo X, Zhao Y, Yuan X. miR-663a inhibits hepatocellular carcinoma cell proliferation and invasion by targeting HMGA2. *Biomed Pharmacother* 2016 Jul;81:431-8.
 49. Zhu Y, Wei W, Ye T, Liu Z, Liu L, Luo Y, Zhang L, Gao C, Wang N, Yu L. Small Molecule TH-39 Potentially Targets Hec1/Nek2 Interaction and Exhibits Antitumor Efficacy in K562 Cells via G0/G1 Cell Cycle Arrest and Apoptosis Induction. *CELL PHYSIOL BIOCHEM* 2016;40(1-2):297-308.
 50. Zhao M, Su Z, Zhang S, Zhuang L, Xie Y, Li X. Suppressive Role of MicroRNA-148a in Cell Proliferation and Invasion in Ovarian Cancer Through Targeting Transforming Growth Factor- β -Induced 2. *Oncol Res* 2016;24(5):353-360.
 51. Zou L, Chai J, Gao Y, Guan J, Liu Q, Du JJ. Down-regulated PLAC8 promotes hepatocellular carcinoma cell proliferation by enhancing PI3K/Akt/GSK3 β /Wnt/ β -catenin signaling. *Biomed Pharmacother* 2016 Dec;84:139-146.
 52. Wang Y, Zhang B, Liu W, Dai Y, Shi Y, Zeng Q, Wang F. Noninvasive bioluminescence imaging of the dynamics of sanguinarine induced apoptosis via activation of reactive oxygen species. *ONCOTARGET* 2016 Apr 19;7(16):22355-67.
 53. Liu C, Qian P, Yang L, Zhang L, Chen C, He M, Lu Y, Feng W, Li M, Zhang Y, Zhong M, Yu Z, Zhou Z. Pubertal exposure to di-(2-ethylhexyl)-phthalate inhibits G9a-mediated histone methylation during spermatogenesis in mice. *Arch Toxicol* 2016 Apr;90(4):955-69.
 54. Zhang Z, Guo M, Zhang J, Du C, Xing Y. Leptin Regulates Tau Phosphorylation through Wnt Signaling Pathway in PC12 Cells. *NEURO SIGNALS* 2016 Oct 24;24(1):95-101.
 55. Guo W, Yu D, Wang X, Luo C, Chen Y, Lei W, Wang C, Ge Y, Xue W, Tian Q, Gao X, Yao W. Anti-inflammatory effects of interleukin-23 receptor cytokine-binding homology region rebalance T cell distribution in rodent collagen-induced arthritis. *ONCOTARGET* 2016 May 31;7(22):31800-13.
 56. Yuan C, Zhu X, Han Y, Song C, Liu C, Lu S, Zhang M, Yu F, Peng Z, Zhou C. Elevated HOXA1 expression correlates with accelerated tumor cell proliferation and poor prognosis in gastric cancer partly via cyclin D1. *J Exp Clin Canc Res* 2016 Jan 21;35:15.
 57. Yang R, Liu M, Liang H, Guo S, Guo X, Yuan M, Lian H, Yan X, Zhang S, Chen X, Fang F, Guo H, Zhang C. miR-138-5p contributes to cell proliferation and invasion by targeting Survivin in bladder cancer cells. *Mol Cancer* 2016 Dec 15;15(1):82.
 58. Ling X, Zhang G, Chen Q, Yang H, Sun L, Zhou N, Wang Z, Zou P, Wang X, Cui Z, Liu J, Ao L, Cao J. Shorter sperm telomere length in association with exposure to polycyclic aromatic hydrocarbons: Results from the MARHCS cohort study in Chongqing, China and in vivo animal experiments. *Environ Int* 2016 Oct;95:79-85.
 59. Li L, Gu M, You B, Shi S, Shan Y, Bao L, You Y. Long non-coding RNA ROR promotes proliferation, migration and chemoresistance of nasopharyngeal carcinoma. *Cancer Sci* 2016 Sep;107(9):1215-22.
 60. Sun L, Wang D, Li X, Zhang L, Zhang H, Zhang Y5. Extracellular matrix protein ITGBL1 promotes ovarian cancer cell migration and adhesion through Wnt/PCP signaling and FAK/SRC pathway. *Biomed Pharmacother* 2016 Jul;81:145-51.
 61. Li ZW, Wu B, Ye P, Tan ZY, Ji YH. Brain natriuretic peptide suppresses pain induced by Bmk I, a sodium channel-specific modulator, in rats. *J Headache Pain* 2016 Dec;17(1):90.
 62. Wang YJ, Liu JZ, Lv P, Dang Y, Gao JY, Wang Y. Long non-coding RNA CCAT2 promotes gastric cancer proliferation and invasion by regulating the E-cadherin and LATS2. *Am J Cancer Res* 2016 Nov 1;6(11):2651-2660.
 63. Li H, Fan J, Sun L, Liu X, Cheng P, Fan H. Functional regeneration of ligament-bone interface using a triphasic silk-based graft. *Biomaterials* 2016 Nov;106:180-92.
 64. Zhang T, Wu Y, Fang Z, Yan Q, Zhang S, Sun R, Khaliq J, Li Y. Low expression of RBMS3 and SFRP1 are associated with poor prognosis in patients with gastric cancer. *Am J Cancer Res* 2016 Nov 1;6(11):2679-2689.
 65. Sun JY, Wang JD, Wang X, Liu HC, Zhang MM, Liu YC, Zhang CH, Su Y, Shen YY, Guo YW, Shen AJ, Geng MY. Marine-derived chromopeptide A, a novel class I HDAC inhibitor, suppresses human prostate cancer cell proliferation and migration. *Acta Pharmacol Sin* 2017 Jan 23.
 66. Zhou ZR, Yang ZZ, Wang SJ, Zhang L, Luo JR, Feng Y, Yu XL, Chen XX, Guo XM. The Chk1 inhibitor MK-8776 increases the radiosensitivity of human triple-negative breast cancer by inhibiting autophagy. *Acta Pharmacol Sin* 2017 Jan 2.
 67. Li Y, Wang LM, Xu JZ, Tian K, Gu CX, Li ZF. *Gastrodia elata* attenuates inflammatory response by inhibiting the NF- κ B pathway in rheumatoid arthritis fibroblast-like synoviocytes. *Biomed Pharmacother* 2017 Jan;85:177-181.
 68. Tang Z, Li J, Shen Q, Feng J, Liu H, Wang W, Xu L, Shi G, Ye X, Ge M, Zhou X, Ni S. Contribution of upregulated dipeptidyl peptidase 9 (DPP9) in promoting tumorigenicity, metastasis and the prediction of poor prognosis in non-small cell lung cancer (NSCLC). *Int J Cancer* 2017 Apr 1;140(7):1620-1632.
 69. Sun Z, Nie Q, Zhang L, Niu R, Wang J, Wang S. Fluoride reduced the immune privileged function of mouse Sertoli cells via the regulation of Fas/FasL system. *Chemosphere* 2017 Feb;168:318-325.
 70. Zhou X, Sun F, Luo S, Zhao W, Yang T, Zhang G, Gao M, Lu R, Shu Y, Mu W, Zhuang Y, Ding F, Xu C, Lu Y. Let-7a is an antihypertrophic regulator in the heart via targeting calmodulin. *Int J Biol Sci* 2017 Jan 1;13(1):22-31.
 71. Liu H, Ai J, Shen A, Chen Y, Wang X, Peng X, Chen H, Shen Y, Huang M, Ding J, Geng M. c-Myc alteration determines the therapeutic response to FGFR

- Inhibitors. *Clin Cancer Res* 2017 Feb 15;23(4):974-984.
72. Shi Y, Sun X, He X. Overexpression of Aristaless-Like Homeobox-4 Inhibits Proliferation, Invasion, and EMT in Hepatocellular Carcinoma Cells. *Oncol Res* 2017 Jan 2;25(1):11-18.
 73. Zhao P, Guan HT, Dai ZJ, Ma YG, Liu XX, Wang XJ. Knockdown of Tripartite Motif-Containing Protein 37 (TRIM37) Inhibits the Proliferation and Tumorigenesis in Colorectal Cancer Cells. *Oncol Res* 2017 Jan 2;25(1):115-122.
 74. Zhou H, Yang C, Chen H, Li X, Li Y, Fan X. A simple G-quadruplex molecular beacon-based biosensor for highly selective detection of microRNA. *Biosens Bioelectron* 2017 Jan 15;87:552-557.
 75. Liu X, Quan B, Tian Z, Xi H, Jia G, Wang H, Zhang L, Liu R, Ma C, Han F, Li H, Yuan F. Elevated expression of KLK8 predicts poor prognosis in colorectal cancer. *Biomed Pharmacother* 2017 Jan 28;88:595-602.
 76. Guo Y, Han B, Luo K, Ren Z, Cai L, Sun L. NOX2-ROS-HIF-1 α signaling is critical for the inhibitory effect of oleanolic acid on rectal cancer cell proliferation. *Biomed Pharmacother* 2017 Jan;85:733-739.
 77. Yang K, Wang Y, Ju Y, Li G, Liu C, Liu J, Liu Q, Wen X, Liu LC. p75 neurotrophin receptor regulates differential mineralization of rat ectomesenchymal stem cells. *Cell Proliferat* 2017 Feb;50(1).
 78. He ZJ, Zhu FY, Li SS, Zhong L, Tan HY, Wang K. Inhibiting ROS-NF- κ B-dependent autophagy enhanced brazilin-induced apoptosis in head and neck squamous cell carcinoma. *Food Chem Toxicol* 2017 Jan 4;101:55-66.
 79. Zheng N, Liu L, Liu WW, Li F, Hayashi T, Tashiro SI, Onodera S, Ikejima T. Crosstalk of ROS/RNS and autophagy in silibinin-induced apoptosis of MCF-7 human breast cancer cells *in vitro*. *Acta Pharmacol Sin* 2017 Feb;38(2):277-289.
 80. Liu J, Li Q, Li R, Ren P, Dong S. MicroRNA-363-3p inhibits papillary thyroid carcinoma progression by targeting PIK3CA. *Am J Cancer Res* 2017 Jan 1;7(1):148-158.
 81. Zhang B, Chang M, Wang J, Liu Y. Neuroglobin functions as a prognostic marker and promotes the tumor growth of glioma via upregulating apoptosis. *Biomed Pharmacother* 2017 Jan 16;88:173-180.
 82. Huang T, Liu R, Fu X, Yao D, Yang M, Liu Q, Lu WW, Wu C, Guan M. Aging Reduces an ERR α -Directed Mitochondrial Glutaminase Expression Suppressing Glutamine Anaplerosis and Osteogenic Differentiation of Mesenchymal Stem Cells. *Stem Cells* 2017 Feb;35(2):411-424.
 83. Mei L, Yuan L, Shi W, Fan S, Tang C, Fan X, Yang W, Qian Y, Hussain M, Wu X. SUMOylation of large tumor suppressor 1 at Lys751 attenuates its kinase activity and tumor-suppressor functions. *Cancer Lett* 2017 Feb 1;386:1-11.
 84. Luo C, Chen S, Xu N, Sai WB, Zhao W, Li YC, Hu XJ, Tian H, Gao XD, Yao WB. Establishment of a fluorescence-based method to evaluate endocytosis of desialylated glycoproteins *in vitro*. *Biomed Pharmacother* 2017 Jan 14;88:87-94.
 85. Lin S, Wang J, Wang L, Wen J, Guo Y, Qiao W, Zhou J, Xu G, Zhi F. Phosphodiesterase-5 inhibition suppresses colonic inflammation-induced tumorigenesis via blocking recruitment of MDSC. *Am J Cancer Res* 2017 Jan 1;7(1):41-52.
 86. Xie Y, Guo H, Wang L, Xu L, Zhang X, Yu L, Liu Q, Li Y, Zhao N, Zhao N, Ye R, Liu X. Human albumin attenuates excessive innate immunity via inhibition of microglial Mincle/Syk signaling in subarachnoid hemorrhage. *Brain Behav Immun* 2017 Feb;60:346-360.
 87. Huang N, Cheng S, Zhang X, Tian Q, Pi J, Tang J, Huang Q, Wang F, Chen J, Xie Z, Xu Z, Chen W, Zheng H, Cheng Y. Efficacy of NGR peptide-modified PEGylated quantum dots for crossing the blood-brain barrier and targeted fluorescence imaging of glioma and tumor vasculature. *Nanomedicine* 2017 Jan;13(1):83-93.
 88. Yin L, Mao Y, Song H, Wang Y, Zhou W, Zhang Z. Vincristine alleviates adriamycin-induced nephropathy through stabilizing actin cytoskeleton. *Cell Biosci* 2017 Jan 3;7:1.
 89. Wang J, Song W, Shen W, Yang X, Sun W, Qu S, Shang R, Ma B, Pu M, Tao K, Dou K, Li H. MicroRNA-200a Suppresses Cell Invasion and Migration by Directly Targeting GAB1 in Hepatocellular Carcinoma. *Oncol Res* 2017 Jan 2;25(1):1-10.
 90. Lan T, Ma W, Hong Z, Wu L, Chen X, Yuan Y. Long non-coding RNA small nucleolar RNA host gene 12 (SNHG12) promotes tumorigenesis and metastasis by targeting miR-199a/b-5p in hepatocellular carcinoma. *J Exp Clin Oncol* 2017 Jan 10;36(1):11.
 91. Huang L, Zhou B, Wu H, Zheng L, Zhao J. Effect of apatite formation of biphasic calcium phosphate ceramic (BCP) on osteoblastogenesis using simulated body fluid (SBF) with or without bovine serum albumin (BSA). *Mat Sci Eng C-Mater* 2017 Jan 1;70(Pt 2):955-961.
 92. Guo J, Hao J, Jiang H, Jin J, Wu H, Jin Z, Li Z. Proteasome activator subunit 3 promotes pancreatic cancer growth via c-Myc-glycolysis signaling axis. *Cancer Lett* 2017 Feb 1;386:161-167.
 93. Li N, Qin S, Xie L, Qin T, Yang Y, Fang W, Chen MH. Elevated Serum Potassium Concentration Alleviates Cerebral Ischemia-Reperfusion Injury via Mitochondrial Preservation. *CELL PHYSIOL BIOCHEM* 2018;48(4):1664-1674.
 94. Qiao JT, Cui C, Qing L, Wang LS, He TY, Yan F, Liu FQ, Shen YH, Hou XG, Chen L. Activation of the STING-IRF3 pathway promotes hepatocyte inflammation, apoptosis and induces metabolic disorders in nonalcoholic fatty liver disease. *Metabolism* 2018 Apr;81:13-24.
 95. Xiang Y, Ye W, Huang C, Yu D, Chen H, Deng T, Zhang F, Lou B, Zhang J, Shi K, Chen B, Zhou M. Brusatol Enhances the Chemotherapy Efficacy of Gemcitabine in Pancreatic Cancer via the Nrf2 Signaling Pathway. *Oxid Med Cell Longev* 2018 Apr 18;2018:2360427.
 96. Wang X, Hu H, Liu H. RNA binding protein Lin28B confers gastric cancer cells stemness via directly binding to NRP-1. *Biomed Pharmacother* 2018 Aug;104:383-389.
 97. Song H, Xu Y, Shi L, Xu T, Fan R, Cao M, Xu W, Song J. LncRNA THOR increases the stemness of gastric cancer cells via enhancing SOX9 mRNA stability. *Biomed Pharmacother* 2018 Dec;108:338-346.
 98. Fang WJ, Wang CJ, He Y, Zhou YL, Peng XD, Liu SK. Resveratrol alleviates diabetic cardiomyopathy in rats by improving mitochondrial function through PGC-1 α deacetylation. *Acta Pharmacol Sin* 2018 Jan;39(1):59-73.
 99. Yao L, Ye Y, Mao H, Lu F, He X, Lu G, Zhang S. MicroRNA-124 regulates the expression of MEKK3 in the inflammatory pathogenesis of Parkinson's disease. *J Neuroinflamm* 2018 Jan 12;15(1):13.
 100. Cao C, Huang Y, Tang Q, Zhang C, Shi L, Zhao J, Hu L, Hu Z, Liu Y, Chen L. Bidirectional juxtacrine EphrinB2/Ephs signaling promotes angiogenesis of ECs and maintains self-renewal of MSCs. *Biomaterials* 2018 Jul;172:1-13.
 101. Wu D, Han R, Deng S, Liu T, Zhang T, Xie H, Xu Y. Protective Effects of Flagellin A N/C Against Radiation-Induced NLR Pyrin Domain Containing 3 Inflammasome-Dependent Pyroptosis in Intestinal Cells. *Int J Radiat Oncol* 2018 May 1;101(1):107-117.
 102. Yang Q, Sun G, Yin H, Li H, Cao Z, Wang J, Zhou M, Wang H, Li J. PINK1 Protects Auditory Hair Cells and Spiral Ganglion Neurons from Cisplatin-induced Ototoxicity via Inducing Autophagy and Inhibiting JNK Signaling Pathway. *Free Radical Bio Med* 2018 May 20;120:342-355.
 103. Luo J, Xiang Y, Xu X, Fang D, Li D, Ni F, Zhu X, Chen B, Zhou M. High Glucose-Induced ROS Production Stimulates Proliferation of Pancreatic Cancer via Inactivating the JNK Pathway. *Oxid Med Cell Longev* 2018 Nov 21;2018:6917206.
 104. Yu B, Liu D, Zhang H, Xie D, Nie W, Shi K, Yang P. Anti-hypertrophy effect of atorvastatin on myocardium depends on AMPK activation-induced miR-143-3p suppression via Foxo1. *Biomed Pharmacother* 2018 Oct;106:1390-1395.
 105. Ma TJ, Zhang ZW, Lu YL, Zhang YY, Tao DC, Liu YQ, Ma YX. CLOCK and BMAL1 stabilize and activate RHOA to promote F-actin formation in cancer cells. *Exp Mol Med* 2018 Oct 4;50(10):130.
 106. Li Y, Jiang X, Zhang Y, Gao Z, Liu Y, Hu J, Hu X, Li L, Shi J, Gao N. Nuclear accumulation of UBC9 contributes to SUMOylation of lamin A/C and nucleophagy in response to DNA damage. *J EXP CLIN ONCOL RES* 2019 Feb 11;38(1):67.
 107. Li F, Tan W, Chen H, Zhou J, Xu M, Yuan G. Up- and downregulation of mature miR-1587 function by modulating its G-quadruplex structure and using small molecules. *Int J Biol Macromol* 2019 Jan 12;127-134.

108. Zhang M, Sun H, Deng Y, Su M, Wei S, Wang P, Yu L, Liu J, Guo J, Wang X, Han X, He Q, Shen L COPI-Mediated Nuclear Translocation of EGFRvIII Promotes STAT3 Phosphorylation and PKM2 Nuclear Localization. *Int J Biol Sci* 2019 Jan 1 15(1):114-126.
109. Zhou X, Chu X, Yuan H, Qiu J, Zhao C, Xin D, Li T, Ma W, Wang H, Wang Z, Wang D Mesenchymal stem cell derived EVs mediate neuroprotection after spinal cord injury in rats via the microRNA-21-5p/FasL gene axis. *Biomed Pharmacother* 2019 Jul 115:108818.
110. Sun K, Tang S, Hou Y, Xi L, Chen Y, Yin J, Peng M, Zhao M, Cui X, Liu M Oxidized ATM-mediated glycolysis enhancement in breast cancer-associated fibroblasts contributes to tumor invasion through lactate as metabolic coupling. *EBioMedicine* 2019 Mar 41:370-383.
111. Huang X, Han X, Huang Z, Yu M, Zhang Y, Fan Y, Xu B, Zhou K, Song L, Wang X, Lu C, Xia Y Maternal pentachlorophenol exposure induces developmental toxicity mediated by autophagy on pregnancy mice. *ECOTOX ENVIRON SAFE* 2019 Mar 169:829-836.
112. Niu Y, Lin Z, Wan A, Chen H, Liang H, Sun L, Wang Y, Li X, Xiong XF, Wei B, Wu X, Wan G RNA N6-methyladenosine demethylase FTO promotes breast tumor progression through inhibiting BNIP3. *Mol Cancer* 2019 Mar 28 18(1):46.
113. Ye W, Zheng C, Yu D, Zhang F, Pan R, Ni X, Shi Z, Zhang Z, Xiang Y, Sun H, Shi K, Chen B, Zhang Q, Zhou M Lipoxin A4 Ameliorates Acute Pancreatitis-Associated Acute Lung Injury through the Antioxidative and Anti-Inflammatory Effects of the Nrf2 Pathway. *Oxid Med Cell Longev* 2019 Nov 6 2019:2197017.
114. Sun Y, Jiang C, Hong H, Liu J, Qiu L, Huang Y, Ye L Effects of hypoxia on cardiomyocyte proliferation and association with stage of development. *Biomed Pharmacother* 2019 Oct 118:109391.
115. Zhang JT, Qin H, Man Cheung FK, Su J, Zhang DD, Liu SY, Li XF, Qin J, Lin JT, Jiang BY, Song Dong, Liao RQ, Qiang N, Yang XN, Tu HY, Zhou Q, Yang JJ, Zhang XC, Zhang YN, Wu YL, Zhong WZ Plasma extracellular vesicle microRNAs for pulmonary ground-glass nodules. *J Extracell Vesicles* 2019 Sep 18 8(1):1663666.
116. Xie Q, Wang Z, Zhou H, Yu Z, Huang Y, Sun H, Bi X, Wang Y, Shi W, Gu P, Fan X. The role of miR-135-modified adipose-derived mesenchymal stem cells in bone regeneration. *Biomaterials* 2016 Jan;75:279-94.
117. Wang X, Hu H, Liu H RNA binding protein Lin28B confers gastric cancer cells stemness via directly binding to NRP-1. *Biomed Pharmacother* 2018 Aug;104:383-389.
118. Song K, Yuan X, Li G, Ma M, Sun J. Long noncoding RNA CASC11 promotes osteosarcoma metastasis by suppressing degradation of snail mRNA. *Am J Cancer Res* 2019 Feb 1 9(2):300-311. eCollection 2019. Retraction in: *Am J Cancer Res*. 2019 Oct 01;9(10):2279.
119. Chang Liu, Jian Wang, Li Zhao, Hui He, Pan Zhao, Zheng Peng, Feiyuan Liu, Juan Chen, Weiqing Wu, Guangsuo Wang, Fajin Dong Knockdown of Thymidine Kinase 1 Suppresses Cell Proliferation, Invasion, Migration, and Epithelial-Mesenchymal Transition in Thyroid Carcinoma Cells *Front Oncol* 2020 Jan 29;9:1475.
120. Guojun Yang, Yixin Yang, Hong Tang, Kai Yang Loss of the clock gene *Per1* promotes oral squamous cell carcinoma progression via the AKT/mTOR pathway *Cancer Sci* 2020 May;111(5):1542-1554.
121. Xiaoying He, Zhaomin Lin, Junya Ning, Na Li, Xiaoling Cui, Baoxiang Zhao, Fanzhen Hong, Junying Miao Promoting TTC4 and HSP70 interaction and translocation of annexin A7 to lysosome inhibits apoptosis in vascular endothelial cells *FASEB J* 2020 Sep;34(9):12932-12945.
122. Qinghai Meng, Yu Li, Tingting Ji, Ying Chao, Jun Li, Yu Fu, Suyun Wang, Qi Chen, Wen Chen, Fuhua Huang, Youran Wang, Qichun Zhang, Xiaoliang Wang, Huimin Bian Estrogen prevent atherosclerosis by attenuating endothelial cell pyroptosis via activation of estrogen receptor α -mediated autophagy *J Adv Res* 2020 Aug 24;28:149-164.
123. Wen-Jin Chen, Li-Hua Chen, Ji Wang, Zhao-Tao Wang, Cui-Ying Wu, Kai Sun, Bo-Yun Ding, Ning Liu, Ru-Xiang Xu LHPP impedes energy metabolism by inducing ubiquitin-mediated degradation of PKM2 in glioblastoma *Am J Cancer Res* 2021 Apr 15;11(4):1369-1390.
124. Dekun Gao, Hui Wu, Huiqun Jie, Min Liang, Dongzhen Yu, Yanmei Feng, Karthikeyan Balasubramanian, Guiliang Zheng, Jun Yang, Jingchun He XIAP inhibits gentamicin-induced hair cell damage and ototoxicity through the caspase-3/9 pathway *Biochem Pharmacol* 2021 Apr;186:114513.
125. Yanyi Li, Qifan Hu, Weishuo Li, Shijing Liu, Kaiming Li, Xiaoyu Li, Junjie Du, Zexuan Yu, Cong Wang, Can Zhang Simultaneous blockage of contextual TGF- β by cyto-pharmaceuticals to suppress breast cancer metastasis *J Control Release* 2021 Aug 10;336:40-53.
126. Licheng Yin, Mengyuan Lv, Xingyang Qiu, Xinyan Wang, Anying Zhang, Kun Yang, Hong Zhou IFN- γ Manipulates NOD1-Mediated Interaction of Autophagy and Edwardsiella piscicida to Augment Intracellular Clearance in Fish *J Immunol* 2021 Aug 15;207(4):1087-1098.
127. Chao Zhou, Mengyu Liu, Xiang Mei, Qian Li, Wenjuan Zhang, Ping Deng, Zhixin He, Yu Xi, Tong Tong, Hui Feng Pi, Yonghui Lu, Chunhai Chen, Lei Zhang, Zhengping Yu, Zhou Zhou, Mindi He Histone hypoacetylation contributes to neurotoxicity induced by chronic nickel exposure in vivo and in vitro *Sci Total Environ* 2021 Aug 20;783:147014.
128. Wenjian Yao, Jianjun Wang, Fanruo Meng, Zibo Zhu, Xiangbo Jia, Lei Xu, Quan Zhang, Li Wei Circular RNA CircPVT1 Inhibits 5-Fluorouracil Chemosensitivity by Regulating Ferroptosis Through MiR-30a-5p/FZD3 Axis in Esophageal Cancer Cells *Front Oncol* 2021 Dec 13;11:780938.
129. Xiaoting Liao, Weikang Zhang, Huijun Dai, Ren Jing, Mengling Ye, Wanyun Ge, Shenglin Pei, Linghui Pan Neutrophil-Derived IL-17 Promotes Ventilator-Induced Lung Injury via p38 MAPK/MCP-1 Pathway Activation *Front Immunol* 2021 Dec 15;12:768813.
130. Lili Zhang, Shilun Li, Juan Li, Yukun Li LncRNA ORLNC1 Promotes Bone Marrow Mesenchymal Stem Cell Pyroptosis Induced by Advanced Glycation End Production by Targeting miR-200b-3p/Foxo3 Pathway *Stem Cell Rev Rep* 2021 Dec;17(6):2262-2275.
131. Yongning Lu, Miao Liu, Nicholas J Tursi, Bin Yan, Xiang Cao, Qi Che, Nianqin Yang, Xi Dong Uropathogenic Escherichia coli Infection Compromises the Blood-Testis Barrier by Disturbing mTORC1-mTORC2 Balance *Front Immunol* 2021 Feb 19;12:582858.
132. Danqing Xin, Tingting Li, Xili Chu, Hongfei Ke, Dexiang Liu, Zhen Wang MSCs-extracellular vesicles attenuated neuroinflammation, synapse damage and microglial phagocytosis after hypoxia-ischemia injury by preventing osteopontin expression *Pharmacol Res* 2021 Feb;164:105322.
133. Xiaotao Xu, Yang Su, Kaixuan Wu, Fan Pan, Aizhong Wang DOCK2 contributes to endotoxemia-induced acute lung injury in mice by activating proinflammatory macrophages *Biochem Pharmacol* 2021 Feb;184:114399.
134. Qing Zhang, Ting Chen, Yun Zhang, Lingxia Lyu, Bohuan Zhang, Chengchen Huang, Xuhao Zhou, Yutao Wu, Zhoubin Li MiR-30c-5p regulates adventitial progenitor cells differentiation to vascular smooth muscle cells through targeting OPG *Stem Cell Res Ther* 2021 Jan 19;12(1):67.
135. Bo Fu, Wei Liu, Cui Zhu, Peng Li, Li Wang, Li Pan, Ke Li, Peiying Cai, Min Meng, Yiting Wang, Anqi Zhang, Wenqiang Tang, Meng An Circular RNA circBCBM1 promotes breast cancer brain metastasis by modulating miR-125a/BRD4 axis *Int J Biol Sci* 2021 Jul 22;17(12):3104-3117.
136. Jin-Zhuo Ning, Kai-Xiang He, Fan Cheng, Wei Li, Wei-Min Yu, Hao-Yong Li, Ting Rao, Yuan Ruan Long Non-coding RNA MEG3 Promotes Pyroptosis in Testicular Ischemia-Reperfusion Injury by Targeting MiR-29a to Modulate PTEN Expression *Front Cell Dev Biol* 2021 Jun 18;9:671613.
137. Jiahui Jin, Qingjian Ou, Zhe Wang, Haibin Tian, Jing-Ying Xu, Furong Gao, Shuqin Hu, Jie Chen, Juan Wang, Jieping Zhang, Lixia Lu, Caixia Jin, Guo-Tong Xu, Jingjun Zhao BMSC-derived extracellular vesicles intervened the pathogenic changes of scleroderma in mice through miRNAs *Stem Cell Res Ther* 2021 Jun 5;12(1):327.
138. Zhanfeng Liang, Qian Zhang, Xue Dong, Zhaoqi Zhang, Hongxia Wang, Jiayu Zhang, Yong Zhao mTORC2 negatively controls the maturation process of medullary thymic epithelial cells by inhibiting the LTBR/RANK-NF- κ B axis *J Cell*

- Physiol 2021 Jun;236(6):4725-4737.
139. Cang Li, Ming-Chong Yang, Pan-Pan Hong, Xiao-Fan Zhao, Jin-Xing Wang Metabolomic Profiles in the Intestine of Shrimp Infected by White Spot Syndrome Virus and Antiviral Function of the Metabolite Linoleic Acid in Shrimp J Immunol 2021 May 1;206(9):2075-2087.
 140. Ying Huang, Shan-He Yu, Wen-Xuan Zhen, Tao Cheng, Dan Wang, Jie-Bo Lin, Yu-Han Wu, Yi-Fan Wang, Yi Chen, Li-Ping Shu, Yi Wang, Xiao-Jian Sun, Yi Zhou, Fan Yang, Chih-Hung Hsu, Peng-Fei Xu Tanshinone I, a new EZH2 inhibitor restricts normal and malignant hematopoiesis through upregulation of MMP9 and ABCG2 Theranostics 2021 May 8;11(14):6891-6904.
 141. Qingbiao Xu, Mingyang Hu, Min Li, Jinxiu Hou, Xianghua Zhang, Ya Gao, Bahram Chachar, Xiang Li Dietary Bioactive Peptide Alanyl-Glutamine Attenuates Dextran Sodium Sulfate-Induced Colitis by Modulating Gut Microbiota Oxid Med Cell Longev 2021 May 8;2021:5543003.
 142. Jin-Jin Zhu, Bin-Yuan Yu, Xiao-Kai Huang, Min-Zhi He, Bin-Wen Chen, Ting-Ting Chen, Huang-Yi Fang, Shang-Qin Chen, Xiao-Qin Fu, Pei-Jun Li, Zhen-Lang Lin, Jiang-Hu Zhu Neferine Protects against Hypoxic-Ischemic Brain Damage in Neonatal Rats by Suppressing NLRP3-Mediated Inflammasome Activation Oxid Med Cell Longev 2021 May 8;2021:6654954.
 143. Jie Chen, Xiao Luo, Mao Liu, Lihui Peng, Zixia Zhao, Chengsong He, Yue He Silencing long non-coding RNA NEAT1 attenuates rheumatoid arthritis via the MAPK/ERK signalling pathway by downregulating microRNA-129 and microRNA-204 RNA Biol 2021 May;18(5):657-668.
 144. Ran Tian, Xin Guan, Hao Qian, Liang Wang, Zhujun Shen, Ligang Fang, Zhenyu Liu Restoration of NRF2 attenuates myocardial ischemia reperfusion injury through mediating microRNA-29a-3p/CCNT2 axis Biofactors 2021 May;47(3):414-426.
 145. Wan Sun, Changchang Hu, Tianyu Wang, Juan Wang, Jieping Zhang, Furong Gao, Qingjian Ou, Haibin Tian, Caixia Jin, Jingying Xu, Jingfa Zhang, Guo-Tong Xu, Lixia Lu Glia Maturation Factor Beta as a Novel Biomarker and Therapeutic Target for Hepatocellular Carcinoma Front Oncol 2021 Nov 2;11:744331.
 146. Pengli Liu, Yilong Cui, Menglin Liu, Bonan Xiao, Jian Zhang, Wanyue Huang, Xuliang Zhang, Miao Song, Yanfei Li Protective effect of mitophagy against aluminum-induced MC3T3-E1 cells dysfunction Chemosphere 2021 Nov;282:131086.
 147. Lingming Kong, Jie Deng, Xiang Zhou, Binbin Cai, Baofu Zhang, Xiaohu Chen, Zongjing Chen, Weiming Wang Sitagliptin activates the p62-Keap1-Nrf2 signalling pathway to alleviate oxidative stress and excessive autophagy in severe acute pancreatitis-related acute lung injury Cell Death Dis 2021 Oct 11;12(10):928.
 148. Jiapeng Bao, Changjian Lin, Xing Zhou, Diana Ma, Lujie Ge, Kai Xu, Safwat Adel Abdo Moqbel, Yuzhe He, Chiyuan Ma, Jisheng Ran, Lidong Wu circFAM160A2 Promotes Mitochondrial Stabilization and Apoptosis Reduction in Osteoarthritis Chondrocytes by Targeting miR-505-3p and SIRT3 Oxid Med Cell Longev 2021 Oct 4;2021:5712280.
 149. Yilu Xu, Lancheng Wei, Shen Tang, Qianqian Shi, Bin Wu, Xiaobo Yang, Yunfeng Zou, Xinhang Wang, Qingqing Ao, Ling Meng, Xuejing Wei, Ning Zhang, Yunqing Li, Chunhua Lan, Muting Chen, Xiyi Li, Cailing Lu Regulation PP2Ac methylation ameliorating autophagy dysfunction caused by Mn is associated with mTORC1/ULK1 pathway Food Chem Toxicol 2021 Oct;156:112441.
 150. Yanning Li, Yitong He, Haoliang Fan, Zhuo Wang, Jian Huang, Gehua Wen, Xiaohan Wang, Qiqian Xie, Pingming Qiu Brain-derived neurotrophic factor upregulates synaptic GluA1 in the amygdala to promote depression in response to psychological stress Biochem Pharmacol 2021 Oct;192:114740.
 151. Pu Li, Yuezong Bai, Boer Shan, Wei Zhang, Zhanjie Liu, Yingjie Zhu, Xiaoya Xu, Qian Chen, Xiujie Sheng, Xiaoyang Deng, Zhengchen Guo, Dadong Zhang, Huaying Wang, Yanan Zhang, Yuanjing Hu Exploration of Potential Diagnostic Value of Protein Content in Serum Small Extracellular Vesicles for Early-Stage Epithelial Ovarian Carcinoma Front Oncol 2021 Sep 15;11:707658.
 152. Jie Huang, Lu Fang, Shenli Zhang, Ying Zhang, Kunlin Ou, Chonggang Wang Long-term exposure to environmental levels of phenanthrene disrupts spermatogenesis in male mice Environ Pollut 2021 Sep 15;285:117488.
 153. Chao Zhang, Xiao-Chang Wu, Sha Li, Lian-Jie Dou, Li Zhou, Feng-Hui Wang, Kai Ma, Dan Huang, Ying Pan, Ji-Jun Gu, Ji-Yu Cao, Hua Wang, Jia-Hu Hao Perinatal low-dose bisphenol AF exposure impairs synaptic plasticity and cognitive function of adult offspring in a sex-dependent manner Sci Total Environ 2021 Sep 20;788:147918.
 154. Yike Wang, Lifeng Dong, Fang Wan, Fangfang Chen, Dianlei Liu, Deqin Chen, Jingpei Long MiR-9-3p regulates the biological functions and drug resistance of gemcitabine-treated breast cancer cells and affects tumor growth through targeting MTDH Cell Death Dis 2021 Sep 22;12(10):861.
 155. Xiaoyi Zhang, Jing Tao, Jia Yu, Ning Hu, Xuanzhe Zhang, Guirong Wang, Jiarui Feng, Xingcheng Xiong, Man Li, Dongqi Chai, Hanjun Li, Yuping Rong, Zhigang Tang, Weixing Wang, Zhiyong Peng, Qiao Shi Inhibition of Notch activity promotes pancreatic cytokeratin 5-positive cell differentiation to beta cells and improves glucose homeostasis following acute pancreatitis Cell Death Dis 2021 Sep 23;12(10):867.
 156. Menglin Liu, Xia Wu, Yilong Cui, Pengli Liu, Bonan Xiao, Xuliang Zhang, Jian Zhang, Zhuo Sun, Miao Song, Bing Shao, Yanfei Li Mitophagy and apoptosis mediated by ROS participate in AICl3-induced MC3T3-E1 cell dysfunction Food Chem Toxicol 2021 Sep;155:112388.
 157. Cheng-Cheng Wang, Lei Du, Hao-Hao Shi, Lin Ding, Teruyoshi Yanagita, Chang-Hu Xue, Yu-Ming Wang, Tian-Tian Zhang Dietary EPA-Enriched Phospholipids Alleviate Chronic Stress and LPS-Induced Depression- and Anxiety-Like Behavior by Regulating Immunity and Neuroinflammation Mol Nutr Food Res 2021 Sep;65(17):e2100009.
 158. Baiwen Chen, Jijia Hu, Xianting Hu, Huifang Chen, Rujuan Bao, Yatao Zhou, Youqiong Ye, Meixiao Zhan, Wei Cai, Huabin Li, Hua-Bing Li DENR controls JAK2 translation to induce PD-L1 expression for tumor immune evasion Nat Commun 2022 Apr 19;13(1):2059.
 159. Yan Zou, Xinhong Sun, Qingshan Yang, Meng Zheng, Olga Shimoni, Weimin Ruan, Yibin Wang, Dongya Zhang, Jinlong Yin, Xiangang Huang, Wei Tao, Jong Bae Park, Xing-Jie Liang, Kam W Leong, Bingyang Shi Blood-brain barrier-penetrating single CRISPR-Cas9 nanocapsules for effective and safe glioblastoma gene therapy Sci Adv 2022 Apr 22;8(16):eabm8011.
 160. Nan Chen, Yong-Jun Liu, Yong-Liang Fan, Xiao-Jin Pei, Yang Yang, Ming-Tao Liao, Jiru Zhong, Na Li, Tong-Xian Liu, Guirong Wang, Yufeng Pan, Coby Schal, Sheng Li A single gene integrates sex and hormone regulators into sexual attractiveness Nat Ecol Evol 2022 Aug;6(8):1180-1190.
 161. Yufan Zheng, Wenrui Sun, Cong Shan, Borui Li, Jiaying Liu, Hao Xing, Qingling Xu, Baiping Cui, Wenjia Zhu, Jia Chen, Liyan Liu, Tian Yang, Ning Sun, Xiaobo Li β -hydroxybutyrate inhibits ferroptosis-mediated pancreatic damage in acute liver failure through the increase of H3K9hbh Cell Rep 2022 Dec 20;41(12):111847.
 162. Jialiang Wang, Kuirong Mao, Xiuxiu Cong, Huizhu Tan, Chenxi Wu, Zheng Hu, Yong-Guang Yang, Tianmeng Sun Nanoparticle delivery of CD40 siRNA suppresses alloimmune responses by inhibiting activation and differentiation of DCs and macrophages Sci Adv 2022 Dec 21;8(51):eabq3699.
 163. Wei Li, Qi Long, Hao Wu, Yanshuang Zhou, Lifan Duan, Hao Yuan, Yingzhe Ding, Yile Huang, Yi Wu, Jinyu Huang, Delong Liu, Baodan Chen, Jian Zhang, Juntao Qi, Shiwei Du, Linpeng Li, Yang Liu, Zifeng Ruan, Zihuang Liu, Zichao Liu, Yifan Zhao, Jianghuan Lu, Junwei Wang, Wai-Yee Chan, Xingguo Liu Nuclear localization of mitochondrial TCA cycle enzymes modulates pluripotency via histone acetylation Nat Commun 2022 Dec 2;13(1):7414.
 164. Anjun Jiao, Chenming Sun, Xin Wang, Lei Lei, Haiyan Liu, Wenhui Li, Xiaofeng Yang, Huiqiang Zheng, Renyi Ding, Kun Zhu, Yanhong Su, Cangang Zhang, Lianjun Zhang, Baojun Zhang DEXD/H-box helicase 9 intrinsically controls CD8+ T cell-mediated antiviral response through noncanonical mechanisms Sci Adv 2022 Feb 11;8(6):eabk2691.
 165. Jie Wang, Junshang Ge, Yan Wang, Fang Xiong, Jiayue Guo, Xianjie Jiang, Lishen Zhang, Xiangying Deng, Zhaojian Gong, Shanshan Zhang, Qijia Yan, Yi He, Xiayu Li, Lei Shi, Can Guo, Fuyan Wang, Zheng Li, Ming Zhou, Bo Xiang, Yong Li, Wei

- Xiong, Zhaoyang Zeng EBV miRNAs BART11 and BART17-3p promote immune escape through the enhancer-mediated transcription of PD-L1 *Nat Commun* 2022 Feb 14;13(1):866.
166. Yi Li, Yu-Chen Li, Xiao-Tian Liu, Lu Zhang, Yi-Hua Chen, Qiong Zhao, Wen Gao, Baolin Liu, Hua Yang, Ping Li Blockage of citrate export prevents TCA cycle fragmentation via Irg1 inactivation *Cell Rep* 2022 Feb 15;38(7):110391.
167. Xiaojie Ma, Yunkun Lu, Ziyu Zhou, Qin Li, Xi Chen, Weiyun Wang, Yan Jin, Zhensheng Hu, Guo Chen, Qian Deng, Weina Shang, Hao Wang, Hongxing Fu, Xiangwei He, Xin-Hua Feng, Saiyong Zhu Human expandable pancreatic progenitor-derived β cells ameliorate diabetes *Sci Adv* 2022 Feb 25;8(8):eabk1826.
168. Jingjing Zhang, Jinxin Qiu, Wenyong Zhou, Jianping Cao, Xuefei Hu, Wenli Mi, Bing Su, Bin He, Ju Qiu, Lei Shen Neuropilin-1 mediates lung tissue-specific control of ILC2 function in type 2 immunity *Nat Immunol* 2022 Feb;23(2):237-250.
169. Zhi Qi, Zihan Xu, Liuzhen Zhang, Yongkang Zou, Jinping Li, Wenyu Yan, Cheng Li, Ningshu Liu, Hong Wu Overcoming resistance to immune checkpoint therapy in PTEN-null prostate cancer by intermittent anti-PI3Ka/ β / δ treatment *Nat Commun* 2022 Jan 10;13(1):182.
170. Jiliang Xia, Jingyu Zhang, Xuan Wu, Wanqing Du, Yinghong Zhu, Xing Liu, Zhenhao Liu, Bin Meng, Jiaojiao Guo, Qin Yang, Yihui Wang, Qinglin Wang, Xiangling Feng, Guoxiang Xie, Yi Shen, Yanjuan He, Juanjuan Xiang, Minghua Wu, Gang An, Lugui Qiu, Wei Jia, Wen Zhou Blocking glycine utilization inhibits multiple myeloma progression by disrupting glutathione balance *Nat Commun* 2022 Jul 11;13(1):4007.
171. Mingyang Li, Xinyang Jin, Tao Liu, Feng Fan, Feng Gao, Shuang Chai, Lihua Yang Nanoparticle elasticity affects systemic circulation lifetime by modulating adsorption of apolipoprotein A-I in corona formation *Nat Commun* 2022 Jul 16;13(1):4137.
172. Jiaye Liu, Yang Wang, Chunyang Mu, Meng Li, Kewei Li, Shan Li, Wenshuang Wu, Lingyao Du, Xiaoyun Zhang, Chuan Li, Wei Peng, Junyi Shen, Yang Liu, Dujiang Yang, Kaixiang Zhang, Qingyang Ning, Xiaoying Fu, Yu Zeng, Yinyun Ni, Zongguang Zhou, Yi Liu, Yiguo Hu, Xiaofeng Zheng, Tianfu Wen, Zhihui Li, Yong Liu Pancreatic tumor eradication via selective Pin1 inhibition in cancer-associated fibroblasts and T lymphocytes engagement *Nat Commun* 2022 Jul 25;13(1):4308.
173. Xueguang Zhang, Lingbo Wang, Yongyi Ma, Yan Wang, Hongqian Liu, Mohan Liu, Lang Qin, Jinghong Li, Chuan Jiang, Xiaojian Zhang, Xudong Shan, Yuliang Liu, Jinsong Li, Yaqian Li, Rui Zheng, Yongkang Sun, Jianfeng Sun, Xiangyou Leng, Yan Liang, Feng Zhang, Xiaohui Jiang, Yihong Yang, Ying Shen CEP128 is involved in spermatogenesis in humans and mice *Nat Commun* 2022 Mar 16;13(1):1395.
174. Cai Han, Lin-Yu Sun, Xue-Qun Luo, Qi Pan, Yu-Meng Sun, Zhan-Cheng Zeng, Tian-Qi Chen, Wei Huang, Ke Fang, Wen-Tao Wang, Yue-Qin Chen Chromatin-associated orphan snoRNA regulates DNA damage-mediated differentiation via a non-canonical complex *Cell Rep* 2022 Mar 29;38(13):110421.
175. Song Xue, Fan Zhou, Tian Zhao, Huimin Zhao, Xuewei Wang, Long Chen, Jin-Ping Li, Shi-Zhong Luo Phase separation on cell surface facilitates bFGF signal transduction with heparan sulphate *Nat Commun* 2022 Mar 2;13(1):1112.
176. Xiao-Ru Ma, Xudong Zhu, Yujie Xiao, Hui-Min Gu, Shuang-Shuang Zheng, Liang Li, Fan Wang, Zhao-Jun Dong, Di-Xian Wang, Yang Wu, Chenyu Yang, Wenhong Jiang, Ke Yao, Yue Yin, Yang Zhang, Chao Peng, Lixia Gao, Zhuoxian Meng, Zeping Hu, Chong Liu, Li Li, Hou-Zao Chen, Yousheng Shu, Zhenyu Ju, Jing-Wei Zhao Restoring nuclear entry of Sirtuin 2 in oligodendrocyte progenitor cells promotes remyelination during ageing *Nat Commun* 2022 Mar 9;13(1):1225.
177. Su'an Tang, Yumei Cao, Zhaopeng Cai, Xiaoyu Nie, Jianzhao Ruan, Zuoqing Zhou, Guangfeng Ruan, Zhaohua Zhu, Weiyu Han, Changhai Ding The lncRNA PILA promotes NF- κ B signaling in osteoarthritis by stimulating the activity of the protein arginine methyltransferase PRMT1 *Sci Signal* 2022 May 24;15(735):eabm6265.
178. Longfei Ma, Dan Xie, Mengcheng Luo, Xiwen Lin, Hengyu Nie, Jian Chen, Chenxu Gao, Shuguang Duo, Chunsheng Han Identification and characterization of BEND2 as a key regulator of meiosis during mouse spermatogenesis *Sci Adv* 2022 May 27;8(21):eabn1606.
179. Jiang Wang, Jing Zhao, Cong Yan, Cong Xi, Chenglin Wu, Jingxiang Zhao, Fengwei Li, Yanhua Ding, Rui Zhang, Shankang Qi, Xingjun Li, Chao Liu, Wanting Hou, Hong Chen, Yiping Wang, Dalei Wu, Kaixian Chen, Hualiang Jiang, He Huang, Hong Liu Identification and evaluation of a lipid-lowering small compound in preclinical models and in a Phase I trial *Cell Metab* 2022 May 3;34(5):667-680.e6.
180. Xueguang Zhang, Rui Zheng, Chen Liang, Haotian Liu, Xiaozhen Zhang, Yongyi Ma, Mohan Liu, Wei Zhang, Yang Yang, Man Liu, Chuan Jiang, Qingjia Ren, Yan Wang, Suren Chen, Yihong Yang, Ying Shen Loss-of-function mutations in CEP78 cause male infertility in humans and mice *Sci Adv* 2022 Oct 7;8(40):eabn0968.

注：更多使用本产品的文献请参考产品网页。

Version 2024.03.12